

CONTROL

A solid-source mass spectrometer data acquisition program for
use with Hewlett-Packard computers

by
Brian D. Marshall

UNITED STATES GEOLOGICAL SURVEY
OPEN-FILE REPORT 89-160

Revision of Mar 27, 1989

This report is preliminary and has not been reviewed for conformity with U.S. Geological Survey editorial standards and nomenclature. Any use of trade names is for descriptive purposes only, and does not imply endorsement by the U.S. Geological Survey. Although this program has been extensively tested, the U.S. Geological Survey cannot guarantee that it will give accurate results for all applications nor that it will work on all computer systems.

Table of Contents

Introduction	1
Requirements	1
Hardware	1
Software	1
Running a sample	1
Getting started	1
Setting the peak positions	2
Run Parameters	2
Data Acquisition	3
Using the General program option	4
Scanning the magnetic field	4
Appendices	6
Algorithms	6
Source code description	6
Variables	6
Subprograms	8
Source code listings	13
CONTROL	13
Subfile	14
KBDM	30
CABDM	33
RBBDM	37
SRBDM	40
IRBDM	45
References	50

Introduction

CONTROL is a computer program which provides a user-friendly environment for operating a simple solid-source mass spectrometer. It was developed for use with modified NBS-type solid-source mass spectrometers. These mass spectrometers have 6-inch radius flight tubes and were originally manufactured in 1968. CONTROL version 1.1 consists of the following program files: CONTROL, Subfile, CABDM, SRBDM, KBDM, RBBDM, and IRBDM.

Requirements

Hardware

CONTROL runs on a Hewlett-Packard series 200 (or 300) microcomputer with at least 512K of RAM and a dual floppy disk drive. Interfaces must be present for controlling the magnetic field and for controlling and reading a digital voltmeter. Naturally, the software will need to be modified for various hardware configurations. The source code described here requires an HPIB printer with graphics capability set to select code 701, a Solartron 7062 Systems Voltmeter with HPIB interface at select code 716, and a GPIO interface for controlling the magnet set to select code 12.

Software

CONTROL requires H-P BASIC version 3.0 (or later) and the following binary extensions must be loaded prior to running CONTROL: IO, HPIB, GPIO, DISC, MS, MAT, CLOCK, TRANS, GRAPH, GRAPHX.

Running a sample

Getting started

To get CONTROL into memory type "LOAD "CONTROL"" and press the Execute key. Then press the Run key. An opening screen will appear which shows the current time and date, a statement containing the version of CONTROL and a soft key menu at the bottom of the screen. In order to prepare the voltmeter for data acquisition press the soft key labelled Setup DVM. The voltmeter will be set to remote control, the local panel will be disabled, and the initialization will be performed. The voltmeter is then set to a 1 second integration time and started tracking.

The five upper soft keys start various data acquisition programs, four programs each dedicated to a single element (potassium, calcium, rubidium, or strontium) and a fifth program for general data acquisition with no mass discrimination correction or spike subtraction. The Erase Disk option allows the user to initialize a floppy disk placed in the left-hand disk drive. The Shut Down option simply clears the screen, sets the voltmeter to a standby mode and stops the program.

As an example, assume you wish to run a strontium sample. Pressing the Strontium key starts the strontium program which immediately asks for a sample name. Any characters are OK and there is a maximum limit of 35 characters for the name. Press Enter to continue.

At this point, the program presents a soft key menu with the options for type of sample being run, unspiked, spiked, or a spike itself. If the sample is spiked then the user will be presented with choices for spike ratios used to correct the data. Note that these choices must be entered ahead of time into the program itself. If you are analyzing a spike then the program will not correct for mass discrimination.

Setting the peak positions

Once the program progresses through the preliminaries, you will be presented with a screen showing the current peak and magnet setting (e.g. 88 = 3580) and a soft key menu as follows:

Step	UP	Set	Field	
Cycle	DOWN	Scan	3580	Escape

These soft keys allow you to center the peaks using the computer in combination with a manual voltage offset device on the mass spectrometer. The **Step** key switches the magnetic field to the next isotope in sequence, returning to the first isotope upon completion of the sequence. The **UP** and **DOWN** keys change the magnetic field one digital step at a time. The knob will change the magnetic field faster if your keyboard has one. Once the peak of interest is centered, use the **Set** key to store the magnetic field value with the peak. The magnetic field is always displayed below the **Field** key; these keys perform no other purpose. The **Cycle** key steps through the peaks automatically waiting 5 seconds on each before switching. It can only be stopped by pressing the **Hold** key which will become active during the Cycle routine. Cycling the magnetic field before centering the peaks eliminates some of the hysteresis in the magnet, allowing for more accurate centering. It is emphasized that centering of the peaks is critical since the program will not adjust the relative peak positions during the run. The program centers the peaks by finding the largest peak, centering it, and then adjusting all other peaks equally. Of course, you may re-center the peaks manually any time during the run. The **Escape** key exits this segment and presents the user with a screen showing the current run parameters.

Run parameters

The Current Run Parameters screen looks similar to the following:

CURRENT RUN PARAMETERS

Total number of blocks= 30

Number of scans per block= 11

Integration times (seconds):

88= 5 87= 7 86= 7 84= 7 85= 9

Delay times (seconds):

Before 88= 5 Before 87= 6 Before 86= 4 Before 84= 4 Before 85= 6

Print means only.

Correct for 87Rb.

This screen always comes up after exiting the peak adjustment routine. If you answer yes to the question "Do you want to change any of these parameters?" then you will be presented with a series of questions asking for new values. Any yes/no question may be answered with a y or Y or n or N. When changing the integration and/or delay times enter all required values separated by commas. Any question may be skipped by pressing the Enter key without entering any value. If changes are made you will have the opportunity to review them before starting or resuming data acquisition.

After reviewing the current run parameters and answering no to the query, you will be presented with a soft key menu allowing you to start data acquisition, abort this analysis, or escape back to the peak adjustment routine. If you select **START** then the program enters its

main portion. If running a strontium or calcium the program checks to see if a disk with a name containing "DATA" is in the left-hand disk drive. You will be required to provide one if necessary. A data file containing the sample name and date is created on this disk; the name for the file is created from the current date and time as "MMYYDDHH" where MM is a two-letter code for the month and YY, DD, and HH are the year, day, and hour, respectively. As each block is finished the means are stored in this file as a precaution against a hardware or software failure. The use of these files will require another program and is left up to the user. The program then flashes `Printer ON LINE?` for five seconds allowing you to turn it on if it is not already on. Data acquisition will then begin.

Data Acquisition

During background measurements, the screen shows the current measured value and time of measurement. A soft key menu is displayed as follows:

Block	Strontium	Field	Below
1 of 20	Baselines	1815	85

All of these keys are for display purposes only; they will do nothing when pressed. They communicate to the user the current block number, the current field value and which baseline is being measured. During background and interference measurements as well as during peak centering the computer may not be interrupted easily by the user.

When the peak tops are being measured the screen shows a plot of the average peak values and displays the current measured value (baselines are subtracted) and time. The soft key menu looks something like this:

Block	Scan	Strontium	Field	
1 of 20	1 of 11	88	3650	ABORT Block

It is similar to the menu displayed during measurement of the baselines, except that it provides a key (ABORT Block) which will exit the peak measurement routine immediately. If the user presses this key, the program will forget the data for that block (including the baselines) and return to the peak adjustment routine. Data acquisition may be resumed after making any desired adjustments to the beam or to the run parameters.

After a block of data is finished (immediately after the last baseline measurement), the computer will display a soft key menu as follows:

	PAUSE	
FINISH	Change Params	ABORT RUN!

The user will have only a few seconds to respond to this menu; the program will start a new block (assuming that the maximum number of blocks has not been reached) if no key is pressed. The PAUSE key simply pauses the program allowing the user to make minor adjustments (typically focussing) to the hardware which is out of the computer's control. Data acquisition can then be resumed; baselines will not be re-measured. The Change Params key also pauses the program, but unlike the PAUSE key, it returns the user to the peak adjustment routine so that the peak centers can be checked and any run parameters can be changed. When data acquisition is resumed, the baselines are re-measured. For this reason, it is a very good idea to choose Change Params when making hardware adjustments which affect the beam intensity. The FINISH key can be pressed if the user decides that enough data have been acquired; the program will quit normally, allowing final editing of the data. The ABORT RUN! key immediately quits the program without any final data output.

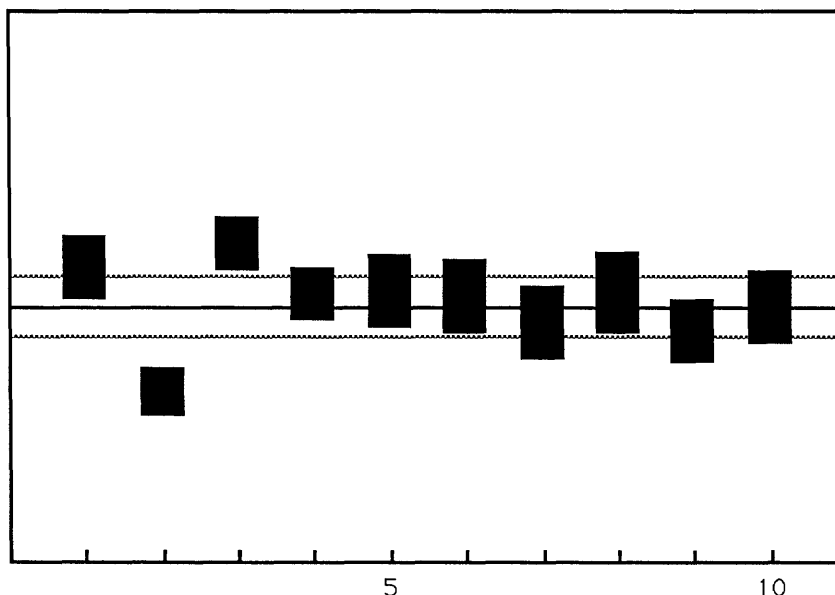


Figure 1. Final plot of an analysis showing the grand mean (solid line) and $\pm 2\sigma$ errors of the mean (dotted lines). Each ratio is plotted as its 2σ error bar.

At the conclusion of an analysis the screen will display a plot (Figure 1) showing the means and 2σ errors for each block. This plot can be dumped to the printer if desired. At this point, the program will ask the user how many blocks to reject. A non-zero answer to this query will be followed by a series of questions asking the user which blocks to reject. The program will continue in this mode allowing the user to re-edit the data over and over (and printing the results) until the user enters "0" for the number of blocks to be rejected.

Using the General program option

The General program allows the user to measure any element but will not perform any mass discrimination correction or spike subtraction; it merely reports raw values. After prompting the user for a sample name, the General program asks the user for an element to be analyzed (version 1.1 supports U, Th, Pb, Re, Rb, K, Sr, and Ca). If the element selected has more than two isotopes the program will ask the user to choose the number of isotopes he/she wishes to measure. The user must then select the isotopes to be measured from a list presented on screen. Note that the reference isotope is the isotope that will be reported in the denominator of the ratios reported. For example, if one wishes to receive data as $^{40}\text{Ca}/^{42}\text{Ca}$ then ^{42}Ca should be chosen as the reference isotope. After choosing the desired isotopes, the General program operates the same as the other options.

Scanning the magnetic field

The Scan soft key is available whenever the user is in the peak adjustment routine. Selecting the Scan key brings up a new set of soft key definitions:

Start Scan	LO field=1450	SCAN UP	
LOG	HI field=3210		Escape

When the scan routine is operating, the voltmeter integration time is set to 0.5 second and the readings are displayed on the screen. The backgrounds should be between 5.2 and 5.3 mV for proper scan display (5 mV is subtracted before plotting). The LO and HI field keys display the lowest and highest magnetic field values to be scanned. The initial values are determined from the current peak settings assuming that the user wishes to scan over the whole mass range being measured. These values may be changed by simply pressing the LO or HI field key and entering a new value. The SCAN UP key indicates that the scan will be from the lowest to the highest field value or up mass. To scan down mass simply press the SCAN UP key, which will then become SCAN DOWN. Note that this key indicates the current setting rather than the setting that will occur if pressed. The LOG key, which operates in a similar manner, indicates that the scan will be plotted with a logarithmic scale. Pressing the LOG key changes it to LINEAR and brings up two more soft key definitions, labelled Min=0V and Max=5V. These two keys refer to the minimum and maximum limits for the plot. They operate just like the LO and HI field keys; note that the values are in volts rather than millivolts.

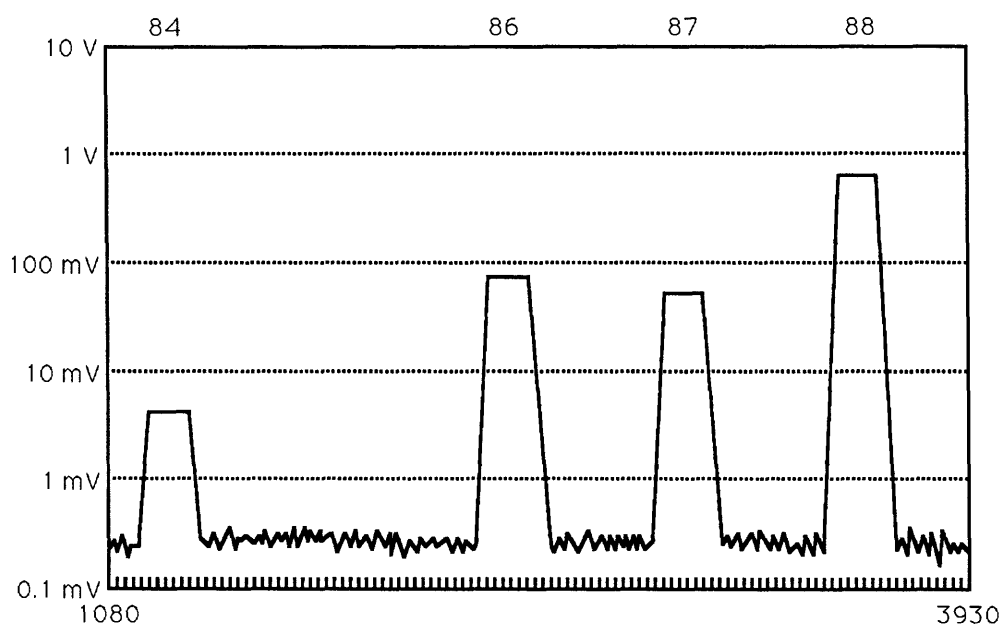


Figure 2. Simulation of plot showing magnetic field scan of a strontium mass spectrum. The mass numbers of the peaks are shown at top and the tick marks at the bottom are shown every 20 digital units.

Once these various options have been selected, pressing the Start Scan key switches the magnet to the first field value (LO or HI field depending on the scan direction) and waits four seconds for the magnet to stabilize before proceeding with the scan. Two soft keys labelled Slower and Faster are active during a scan; they lower or raise the scan speed by factors of two, limited by 1 and 32 as the lowest and highest number of digits per step. The default speed is 4 digits per step. The magnetic field value is displayed in a soft key label during the scan. The Escape key returns the user to the peak adjustment routine. If a scan is completed (the Escape key is not pressed) a soft key labelled Dump Graphics allows the scan to be printed on the printer, (Figure 2).

Appendices

Algorithms

Many users express an interest in learning about the methods used to acquire and reduce the data. This section explains in simple terms the various algorithms used in CONTROL. During data acquisition the backgrounds (measured on each side of each peak) and peak tops are integrated for a given length of time, typically 5 seconds. After every integration the data is edited for outliers. First, data outside 3σ are rejected on the basis of their extremely low probability. Then Chauvenet's criterion is used to reject any data with a probability of occurrence of less than 0.5 (Taylor, 1982). Since backgrounds are measured on both sides of a peak, these two measurements are averaged to obtain the background for a given peak. Backgrounds are only measured before and after each block of peak top measurements, so after a block of data (backgrounds, peak tops, backgrounds) is obtained the backgrounds are extrapolated linearly to the times of measurement of each peak and subtracted from the peak top values to obtain the net peak values. The net peak values are used to obtain ratios using Dodson's (1978) algorithm which corrects for first and second order fluctuations of the ion beam. These raw ratios are then checked for outliers by fitting the ratios to a straight line (in case of varying mass discrimination during a block) and editing them twice as described above for the integrations. The errors estimated for each ratio are used to estimate the internal error for the block. For calcium and strontium measurements, mass discrimination is corrected using an exponential law (Wasserburg *et al.*, 1981). This complicates the spike subtraction routine which must use an iterative calculation to estimate the mass discrimination and sample/spike ratio.

Source code description

Since the program files are not extensively commented, this section describes briefly the meaning of all variables and the operation of all subprograms. These subprograms contain the essential data acquisition, analysis and display routines; they are stored in a separate file called "Subfile" and are loaded into memory by the main programs. The element-specific subroutines such as the default field values, spike ratios, and mass discrimination correction are contained in the program file for that element.

Variables. The following variables are declared at the beginning of each module (KBDM, CABDM, RBBDM, SRBDM, and IRBDM). They are thus global variables; additional variables are allocated by the various subprograms as discussed below.

<u>Name</u>	<u>Type</u>	<u>Usage</u>
A\$	string[4]	="DATA" first 4 characters of disk label of data storage disk
Block\$	string[8]	key label specifying current block and total blocks
Elem\$	string[2]	abbreviated name of element
Element\$	string[10]	full name of element
File\$	string[8]	name for data storage file on disk
N\$	string[6]	disk label of data storage disk
Sample\$	string[35]	sample designation
Traw\$	string array[8]	times of each ratio measurement (HH:MM:SS)
B	integer	loop counter
Block	integer	current block number
Blockmax	integer	total number of blocks to be acquired
Center	integer	peak number of peak to be used for centering

Check	integer(T/F)	true if any run parameters in Parameters sub changed
Delay	integer array	delay times in seconds for each peak
Edit1	integer	number of blocks to be edited
Factor	integer(T/F)	true if total printout desired
False	integer(T/F)	=0; boolean value for false
Halt	integer(T/F)	true if run is paused between blocks
I	integer	loop counter
Index1	integer	subscript of array to be plotted in Summ_plot sub
Index2	integer	subscript of array where deviation is stored
Interference	integer(T/F)	true if an interfering peak is to be corrected for
Intn	integer array	integration times for peaks in seconds
Iter	integer	number of iterations used in spike subtraction routine
Mass	integer array	mass numbers of peaks
Numpeak	integer	number of peaks to be measured
Numsel	integer	Number of isotopes to select from (General option only)
P_mass	integer array	possible mass numbers (General option only)
P_pot	integer array	possible peak field values (General option only)
Params	integer(T/F)	true if run paused and user requested run parameters
Pot	integer array	magnetic field values for peaks
Prog	integer	type of sample (program) to be run (spiked vs unspiked)
Rejected	integer array	number of values rejected from each ratio
Scan	integer	number of scans per block
Sel	integer	number of next isotope to measure (General option only)
Sp	integer	number corresponding to spike selected by user
Timbegin	integer	time at start of data acquisition (minutes)
Timfirst	integer	time at start of block since start of run (minutes)
Timlast	integer	time at finish of block since start of run (minutes)
True	integer(T/F)	=1; boolean value for true
Alpha	real	approximation of spike to sample ratio
Amu	real array	atomic mass of each isotope
Avetime	real array	average time of measurement for 4 peaks used for a ratio
Baseline	real array	values of baselines before and after each block
Beta	real	approximation of normalization ratio
Disc	real	discrimination per mass unit
Ipeak	real array	value of peak used to correct for interference
Itime	real array	time of measurement of interference peak
Mean	real array	block to block average for each ratio
Meancorr	real array	discrimination corrected ratio means within block
Meanraw	real array	raw ratio means within block
Netpeak	real array	baseline corrected peak values
P_exp	real	incremental mass discrimination exponent
Peak	real array	peak values
Percent_dev	real array	percent deviations (1 sigma) for each ratio within block
Qpeak	real array	peak values from quick scan prior to centering peaks
Rcorr	real array	corrected ratios for all blocks
Rraw	real array	raw ratios for each block
Rspike	real array	spike ratios for selected spike
Rstore	real array	array used for storing data on disk
Scantime	real	estimate of time for each block of peak scans in minutes
Sigma	real array	block to block 2-sigma error estimates for each ratio

Sigma_int	real	block to block estimate of total internal error (2-sigma)
Sum_p	real	total mass discrimination exponent
Tcorr	real array	temporary corrected ratios used in spike subtraction routine
Timebase	real array	times of measurement of baselines
Timep	real array	times of measurement of peaks
X_gdu_max	real	horizontal dimension of screen in graphic units
Y_gdu_max	real	vertical dimension of screen in graphic units

Subprograms. The subprograms are listed here in alphabetical order as they are in the Subfile module. The names of subprograms are bold and the variable names are shown in italics. each subprogram is described in three parts: a sentence describing its purpose, another sentence specifying the context in which it is called, and a paragraph describing its operation.

Name: **Center_peaks**
 Purpose: centers the peaks by adjusting their magnetic field values
 Context: called by main program before measuring peaks
 Description: This subprogram first calculates a value (*Side*) by which to adjust the magnetic field so that a position on the side of the peak is obtained. This value is calibrated empirically. It then loops through the peaks in order waiting 4 seconds between each jump. It stores the peak values (*Qpeak*) for use in scaling the plotting routines. This loop also reduces hysteresis effects. During the **Center_peaks** routine the integration time of the voltmeter is changed to 0.5 second. The main loop repeats until an acceptable center position on the centering peak is obtained. This is done by measuring values on the low side, the top, and the high side of the centering peak. A case statement then uses the ratio of the differences from the top to each side to determine how much to adjust the field closer to the center position. These amounts are calibrated empirically. The subprogram then adjusts the magnetic field values for each peak (*Poi*) for the total amount of adjustment calculated for the centering peak. After setting the magnetic field to the first peak position it returns control to the main program.

Name: **Clear_crt**
 Purpose: clears both the alpha and graphics screens
 Context: called by many other contexts

Name: **Dump**
 Purpose: dumps graphics screen to the printer
 Context: called by soft key from certain subprograms

Name: **Edit**
 Purpose: edits blocks and calculates new means and sigmas after a run
 Context: called by main program

Description: The main program, after asking the user how many blocks he/she wishes to edit from the run, calls **Edit** to accomplish the task. **Edit** queries the user as to which blocks he/she wants to edit and then calculates new weighted means. **Edit** then calls **Grandstats** to calculate new external 2-sigmas and the value for the internal 2-sigma for the remaining blocks. **Print_means** is called to print out the results. Note that **Edit** does not change the corrected ratios allowing the user to re-edit the data as many times as necessary, each time beginning with the whole data set.

Name: **Edit_routine**

Purpose: edits each integration for outliers
Context: called by **Measure_peaks**, **Measure_zeros**, and **Measure_inter**
Description: Peak and baseline values are obtained by integrating over a period of time, during which the voltmeter is sending readings to the computer. **Edit_routine** takes these values and returns the average after rejecting outliers. Outliers are rejected at two levels of significance. First, it is assumed that all acceptable values fall within 3 standard deviations of the mean value. Values outside 3 sigma are rejected. Second, the routine checks for values with a probability of occurrence of less than 0.5 (Chauvenet's criterion). In order to do this the program calculates a sigma multiplier (*Sigmult*) such that values outside *Sigmult* standard deviations will have a probability of occurrence of less than 0.5 and therefore can be rejected. *Sigmult* is calculated from a best-fit curve to data from probability tables. **Edit_routine** returns the mean after rejecting all outliers found. If the number of values either before or after rejection falls to 1, **Edit_routine** does not try to reject data but returns the mean.

Name: **Field_values**
Purpose: allows user to set magnetic field values, cycle and scan the magnet field
Context: called by main program
Description: **Field_values** is called at the beginning of a run to get the peak centers from the user. It can also be invoked during a run by aborting a block or by choosing to change the run parameters between blocks. This subprogram can cycle the magnet through the peak settings indefinitely and calls **Magnet_scan** to scan the magnetic field. The computer does not know where the peaks are initially, except that the approximate peak field values are retained as default values from the main program. The user must center the peaks manually using the computer before data acquisition. The program can be in this context indefinitely, during which time the magnetic field can be adjusted, cycled or set using soft keys. The knob will also change the magnetic field. When this subprogram is exited (via a soft key) the magnetic field values should be set for all peaks. Because an interference peak may be too small to easily center manually, its position is calculated based on the settings for the other peaks.

Name: **Final_print**
Purpose: prints the corrected ratios for all blocks at the end of a run
Context: called by the main program
Description: **Final_print** consists of a single loop through the number of blocks of data in order to print out the corrected ratios (*Rcorr*) for each block including the percent error and block number.

Name: **Grandstats**
Purpose: calculates block to block statistics
Context: called by main program and **Edit**
Description: **Grandstats** calculates the block to block or external statistics at the 2-sigma level. Individual blocks are weighted by the inverse square of their estimated percent error and weighted means are calculated. External errors are reported as twice the standard error of the mean.

Name: **Identify**
Purpose: identifies the element to be run and prompts the user for a sample name
Context: called by main program

Name: **Linear_regress**
Purpose: edits each ratio for outliers
Context: called by **Statistics**

Description: **Linear_regress** calculates the means and standard deviations for the raw ratios after rejecting outliers based on Chauvenet's criterion (see **Edit_routine**). It fits the raw ratios to a straight line as a function of their time of measurement because monotonically decreasing or increasing values for a ratio are frequently obtained due to changing mass discrimination during a block. The mean value is the same as would be calculated using a simpler routine, but the error will be more accurate (and lower) than would be calculated otherwise.

Name: **Magnet_scan**

Purpose: scans the magnetic field following users directions

Context: called by **Field_values**

Description: **Magnet_scan** allows the user to scan the magnetic field over a limited gauss range covering the peaks to be measured. The scans are output to the screen and can be dumped to the printer if desired. The user may select (via soft keys) the range of field values to be scanned, whether to scan up field or down field, to draw the scan in log or linear units and if linear then what scale to draw the plot at. The speed of the scan (in field units per integration time) can also be changed during the scan. The integration time of the voltmeter is changed to 0.5 second until the program leaves the **Magnet_scan** routine. The program remains in the **Magnet_scan** routine until the user escapes by pressing a soft key.

Name: **Measure_inter**

Purpose: measures the peak of an interfering element

Context: called by main program

Description: **Measure_inter** measures an isotope peak for an interfering element. It is only implemented for measuring the rubidium interference when measuring strontium samples. The interfering peak is measured after the baselines but before the peak tops. The interfering peak is measured only once between blocks as are the baselines (except for the first block or whenever the user pauses the program to change parameters). For a given block, the interfering peak value and its time of measurement are *Ipeak(1)* and *Itime(1)* for before the block and *Ipeak(2)* and *Itime(2)* after the block. Before each block **New_zeros** is called to set the baselines and interfering peak values to the values measured after the previous block.

Name: **Measure_peaks**

Purpose: measures the peak tops

Context: called by main program

Description: **Measure_peaks** measures the peak tops integrating each for the time specified by the user (stored in *Intn*). Standard graphics statements allow the peak values to be shown on the screen. Approximate baseline subtracted values are shown on the screen and plotted as averages. A soft key is enabled during acquisition of peak tops to allow the user to abort the block. The program then invokes the **Showdefaults** subroutine which shows the current run parameters and calls **Field_values** to allow the user to make any adjustments to the mass spectrometer as necessary.

Name: **Measure_zeros**

Purpose: measures the baselines on both sides of each peak including interfering peak

Context: called by main program

Description: **Measure_zeros** first calculates the magnetic field offset from the peak at which to measure baselines for the peaks. The value in *Del* is the offset in mass

units. On the NBS 6" mass spectrometers this value is about 0.3 mass units for potassium and calcium and nearly 0.5 mass units for rubidium and strontium. The offset in magnetic field units is calculated from the positions (field values) and masses of the first two peaks. Each baseline is integrated for twice the time as its associated peak after a delay equal to the delay before measuring the first peak. To minimize large jumps in the magnetic field, the baseline data is acquired either continuously down mass or up mass depending on the order of peak measurement. If an interfering peak is being measured, its baselines are always measured first.

Name: **New_zeros**

Purpose: sets values of baselines measured after peak tops to values for before peak tops

Context: called by main program

Description: **New_zeros** simply takes the values of the baselines (*Baseline*) and their times of measurement (*Timebase*) from after the last block and sets the values for before the next block equal to them. This allows the baselines to be measured only once between blocks. The interference peak is handled the same way.

Name: **Parameters**

Purpose: displays current run parameters and allows user to change them

Context: called by main program from Showdefaults subroutine

Description: **Parameters** allows the user to view and change the following: total number of blocks to be acquired, the number of scans per block, the peak integration times and delay times, and whether or not all raw ratios are to be printed out or just the means. If strontium is being analyzed, the user can decide whether or not rubidium is to be measured to correct for interference. If potassium or rubidium is being analyzed then the user may change the mass discrimination correction to be applied to the data.

Name: **Plot**

Purpose: sets up the screen to plot the peak tops during a block

Context: called by main program

Description: **Plot** uses standard graphics commands to draw axes and label them for the plot of peak intensities vs time since start of run. The maximum value for the graph is calculated from the maximum peak value obtained by the **Center_peaks** subprogram.

Name: **Print_means**

Purpose: prints block to block means and external errors

Context: called by main program (Printout3 and Summary subroutines) and **Edit**

Description: Once two blocks have been acquired, **Print_means** prints the weighted means and external 2-sigma errors of the means. It is also called after data acquisition to print the final means and is called again to print the means obtained after editing the data.

Name: **Print_out1**

Purpose: prints peak and baseline intensities, magnet field values, and raw ratios

Context: called by main program

Description: **Print_out1** prints the following information after each block: baseline intensities (mV) both before and after the block, initial and final peak intensities (mV), magnet field values, and if the user requested the raw ratios and times of measurement of each ratio.

Name: **Print_out2**

Purpose: prints means of raw ratios, percent deviations and number rejected

Context: called by main program

- Description: **Print_out2** prints the following information after each block: means of the raw ratios, percent deviations (1-sigma) of the means, and the number of ratios that were rejected.
- Name: **Ratios**
- Purpose: calculates the net peak intensities, raw ratios and the times of measurement
- Context: called by main program
- Description: **Ratios** first calculates the rate of change of the baselines for each peak. Using this rate, the baselines are extrapolated linearly to the times of measurement of each peak and subtracted from the peak values to obtain the net peak values (*Netpeak*). Any interfering peak correction is made in a similar way and subtracted from the appropriate *Netpeak* value if significant. The rate of change of the net peaks is then calculated. Using this rate each two pairs of peak measurements is extrapolated linearly to their mean time of measurement (*Avertime*) and their ratio is obtained at this time. This method is used because it inherently corrects for second order rates of change of the peaks (Dodson, 1978). There will be one less ratio obtained than the number of scans measured.
- Name: **Start_up**
- Purpose: initializes variables and starts data acquisition
- Context: called by main program
- Description: **Start_up** asks the user to be sure the printer is on line and then prints the sample name and the date. *Timbegin* is initialized to the current time and *Block* is set to 1.
- Name: **Statistics**
- Purpose: calculates the mean and percent deviations of the raw ratios
- Context: called by main program
- Description: **Statistics** first normalizes the times of measurement by subtracting a constant time from each. Then it calls **Linear_regress** to obtain the regression of the raw ratios as a function of their time of measurement. It then calculates the standard deviation (1-sigma) of the means and returns the percent deviations as well as the mean values.
- Name: **Summ_plot**
- Purpose: plots the corrected ratio of interest for each block on the screen
- Context: called by Summary subroutine of main program
- Description: **Summ_plot** starts by obtaining the corrected ratios for the main ratio of interest and its associated error and storing these as well as the associated mean and error values in temporary variables. It then plots a graph on the screen with an unlabeled y-axis corresponding to the value of the corrected ratio and an x-axis with the block numbers labeled. The minimum value for the plot is calculated from the minimum value of the corrected ratio minus the maximum error. The maximum value for the plot is similarly obtained from the maximum ratio and the maximum error. The mean is plotted as a solid horizontal line and 2-sigma error of the mean is shown by two dashed horizontal lines. Soft keys allow the user to dump the graphic screen to the printer or continue. This plot is very helpful in examining the data during editing.
- Name: **Weird_event**
- Purpose: beeps when certain soft keys are pressed
- Context: called by many other contexts
- Description: **Weird_event** simply returns a beep sound when a soft key is pressed that does not do anything. Since some of the soft key labels are used to

communicate information to the operator, these keys are not supposed to trigger an event. The **Weird_event** subprogram provides a way to effectively disable a soft key but still allow its label to appear on the screen.

Source Code Listings. The complete listings of the source code for CONTROL 1.1 are shown below. These are shown primarily for documentation purposes. The program files are available on disk directly from the author at U.S. Geological Survey, MS 963, Denver, CO 80225.

CONTROL

```

10  REM CONTROL      Version 1.1                Brian D. Marshall  7/88
20  REM This program and all associated modules constitute CONTROL.
30  PRINTER IS CRT
40  DISP CHR$(128);
50  INTEGER True,False
60  DIM Name$(6)
70  False=0
80  True= NOT False
90  GCLEAR
100 OUTPUT KBD;CHR$(255)&CHR$(75);
110 PRINT USING "10(/)"
120 PRINT "USGS Computer Program 'CONTROL'. Version 1.100;"
130 PRINT "Technical Contact: Brian Marshall."
140 ON CYCLE 1,2 GOSUB Update
150 ON KEY 9 LABEL "Shut Down" GOTO Exit
160 ON KEY 0 LABEL "Potassium" GOTO Getk
170 ON KEY 1 LABEL "Calcium" GOTO Getca
180 ON KEY 2 LABEL "Rubidium" GOTO Getrb
190 ON KEY 4 LABEL "General" GOTO Getir
200 ON KEY 3 LABEL "Strontium" GOTO Getsr
210 ON KEY 6 LABEL "Setup DVM" GOSUB Getset
220 ON KEY 5 LABEL "Erase Disk" GOSUB Initdisk
230 WHILE True
240   END WHILE
250 Update:
260   PRINT TABXY(1,1);DATE$(TIMEDATE)
270   PRINT
280   PRINT TIME$(TIMEDATE)
290   RETURN
300 Getk:
310   LOAD "KBDM"
320 Getca:
330   LOAD "CABDM"
340 Getrb:
350   LOAD "RBBDM"
360 Getir:
370   LOAD "IRBDM"
380 Getsr:
390   LOAD "SRBDM"
400 Getset:
410   DISP "Setting up digital voltmeter... please wait."
420   DISABLE
430   REMOTE 716
440   LOCAL LOCKOUT 7
450   OUTPUT 716;"INITIALISE"
460   WAIT 3
470   OUTPUT 716;"LITERALS OFF"
480   OUTPUT 716;"ITIME USER=1"
490   OUTPUT 716;"TRACK ON"
500   OUTPUT 716;"RESUME ON"
510   ENABLE
520   DISP CHR$(12);
530   RETURN

```

```

540 Initdisk:!
550  DISABLE
560  ON ERROR GOTO Done
570  DISP "Insert a disk to be ERASED into LEFT drive and press Continue."
580  PAUSE
590  DISP "Initializing disk...please wait."
600  INITIALIZE ":",700,0"
610  INPUT "Name (6 characters beginning with a letter) for this disk?",Name$
620  PRINT LABEL Name$ TO ":",700,0"
630 Done:!
640  OFF ERROR
650  ENABLE
660  DISP CHR$(12);
670  RETURN
680 Exit:!
690  OFF CYCLE
700  OUTPUT KBD;CHR$(255)&CHR$(75);
710  OFF KEY
720  OUTPUT 716;"TRACK OFF"
730  OUTPUT 716;"DISPLAY [          ]"
740  END

```

Subfile

```

10  REM Subfile                      Brian D. Marshall  4/88
20  SUB Center_peaks(Qpeak(*),INTEGER Numpeak,Interference,Mass(*),Pot(*))
30  OPTION BASE 1
40  INTEGER Delta,J,Sumdelta,K,Side,Center
50  REAL V(3)
60  DIM F$(4)
70  DISP CHR$(128);
80  OUTPUT 716;"ITIME USER=.5"
90  FOR J=1 TO Numpeak+Interference
100  OUTPUT 12 USING "#,W";Pot(J)
110  DISP "Quick scan: ";Mass(J);" = ";Pot(J)
120  WAIT 4
130  OUTPUT 716;"HISTORY CLEAR"
140  ENTER 716;Qpeak(J)
150  NEXT J
160  FOR J=1 TO Numpeak+Interference
170  IF MAX(Qpeak(*))=Qpeak(J) THEN Center=J
180  NEXT J
190  OUTPUT 12 USING "#,W";Pot(Center)
200  DISP CHR$(129);
210  Side=PROUND(ABS(Pot(1)-Pot(2))/ABS(Mass(1)-
Mass(2)))/450*(.684*Mass(Center)+12.6),0)
220  WAIT 4
230  Sumdelta=0
240  REPEAT
250  DISP "Centering mass ";Mass(Center);"      Field= ";Pot(Center);"      New Field=
";Pot(Center)+Sumdelta
260  FOR J=-1 TO 1
270  OUTPUT 12 USING "#,W";Pot(Center)+Sumdelta+J*Side
280  WAIT 3
290  OUTPUT 716;"HISTORY CLEAR"
300  ENTER 716;V(J+2)
310  NEXT J
320  SELECT (V(2)-V(1))/(V(2)-V(3))
330  CASE >4
340  Delta=PROUND(Side/40,0)*4
350  CASE <.25
360  Delta=PROUND(Side/40,0)*(-4)
370  CASE >2
380  Delta=PROUND(Side/40,0)*2
390  CASE <.5
400  Delta=PROUND(Side/40,0)*(-2)

```

```

410     CASE >1.2
420         Delta=PROUND(Side/40,0)
430     CASE <.83
440         Delta=PROUND(Side/40,0)*(-1)
450     CASE ELSE
460         Delta=0
470     END SELECT
480     Sumdelta=Sumdelta+Delta
490     UNTIL Delta=0
500     FOR K=1 TO Numpeak+Interference
510         Pot (K)=Pot (K)+Sumdelta
520     NEXT K
530     OUTPUT 12 USING "#,W";Pot (Center)
540     OUTPUT 716;"ITIME USER=1"
550     DISP CHR$(12);
560 SUBEND
570 SUB Clear_crt
580     OUTPUT KBD;CHR$(255)&CHR$(75);
590     OUTPUT KBD;CHR$(255)&CHR$(84);
600     GCLEAR
610 SUBEND
620 SUB Dump
630     DISABLE
640     DISP CHR$(128);"Dumping graphics...please wait."
650     DUMP GRAPHICS
660     ENABLE
670     PRINT
680     DISP CHR$(12);
690 SUBEND
700 SUB Edit (R(*),Mean(*),Sigma(*),Sigma_int,INTEGER Nedit,Nratio,Blockmax)
710     OPTION BASE 1
720     ALLOCATE W(Blockmax),Rtemp(7,Blockmax)
730     INTEGER N,Bnum,K,Nblock
740     FOR N=1 TO Blockmax
750         W(N)=N
760     NEXT N
770     FOR N=1 TO Nedit
780         INPUT "Number of next block to edit?",Bnum
790         PRINT Bnum;
800         W(Bnum)=0
810     NEXT N
820     PRINT "rejected."
830     Nblock=0
840     FOR N=1 TO Blockmax
850         IF W(N)=0 THEN Goto
860         Nblock=Nblock+1
870         FOR K=1 TO Nratio
880             Rtemp(K,Nblock)=R(K,N)
890         NEXT K
900         Rtemp(7,Nblock)=R(7,N)
910 Goto: !
920     NEXT N
930     CALL Grandstats (Rtemp(*),Mean(*),Sigma(*),Sigma_int,Nratio,Nblock)
940     CALL Print_means (Mean(*),Sigma(*),Sigma_int,Nblock,Nratio)
950 SUBEND
960 SUB Edit_routine (Values(*),Cmean,INTEGER Num,Rej)
970     INTEGER K,J
980     REAL Sigmult,Sumvalues,Sumd,Stdd,D
990     Rej=0
1000     GOSUB Mean_sd
1010     Sigmult=3
1020     GOSUB Edit
1030     GOSUB Mean_sd
1040     Sigmult=1.16+(Num-Rej)*.0988-(Num-Rej)^2*.00214
1050     GOSUB Edit
1060     GOSUB Mean_sd

```

```

1070 SUBEXIT
1080 Mean_sd:
1090 Sumvalues=0
1100 FOR K=1 TO Num-Rej
1110 Sumvalues=Sumvalues+Values(K)
1120 NEXT K
1130 Cmean=Sumvalues/(Num-Rej)
1140 IF Num-Rej<2 THEN SUBEXIT
1150 Sumd=0
1160 FOR K=1 TO Num-Rej
1170 Sumd=Sumd+(Values(K)-Cmean)^2
1180 NEXT K
1190 Stdd=SQR(Sumd/(Num-Rej-1))
1200 RETURN
1210 Edit:
1220 J=0
1230 FOR K=1 TO Num-Rej
1240 D=Values(K)-Cmean
1250 IF ABS(D)>Sigmult*Stdd THEN
1260 Rej=Rej+1
1270 ELSE
1280 J=J+1
1290 Values(J)=Values(K)
1300 END IF
1310 NEXT K
1320 RETURN
1330 SUBEND
1340 SUB Field_values(X_gdu_max,Y_gdu_max,INTEGER Numpeak,Interference,Mass(*),Pot(*))
1350 INTEGER Field,True,I
1360 REAL Dvm
1370 DIM F$(4)
1380 CALL Clear_crt
1390 True=-1
1400 OFF KEY
1410 DISP CHR$(128);
1420 ON KEY 9 LABEL "Escape",2 GOTO Done
1430 ON KEY 0 LABEL "Step",2 GOSUB Step
1440 ON KEY 1 LABEL "UP",2 GOSUB Inc
1450 ON KEY 6 LABEL "DOWN",2 GOSUB Dec
1460 ON KEY 2 LABEL "Set",2 GOSUB Set
1470 ON KEY 5 LABEL "Cycle",2 GOSUB Cycle
1480 ON KEY 3 LABEL "Field",2 CALL Weird_event
1490 ON KEY 7 LABEL "Scan",2 GOSUB Scan
1500 I=1
1510 OUTPUT 12 USING "#,W";Pot(I)
1520 DISP USING "3D,"" = "",4Z";Mass(I);Pot(I)
1530 Field=Pot(I)
1540 GOSUB Update_field
1550 ON KNOB .1,3 GOSUB Turn
1560 WHILE True
1570 END WHILE
1580 Step:
1590 I=I MOD (Numpeak+Interference)+1
1600 OUTPUT 12 USING "#,W";Pot(I)
1610 DISP USING "3D,"" = "",4Z";Mass(I);Pot(I)
1620 Field=Pot(I)
1630 GOSUB Update_field
1640 RETURN
1650 Inc:
1660 Field=Field+1
1670 OUTPUT 12 USING "#,W";Field
1680 GOSUB Update_field
1690 RETURN
1700 Dec:
1710 Field=Field-1
1720 OUTPUT 12 USING "#,W";Field

```

```

1730   GOSUB Update_field
1740   RETURN
1750 Set:~
1760   Pot(I)=Field
1770   DISP USING "3D,"" = "",4Z";Mass(I);Pot(I)
1780   RETURN
1790 Turn:~
1800   Field=Field+KNOBX
1810   IF Field<0 THEN Field=0
1820   IF Field>4095 THEN Field=4095
1830   OUTPUT 12 USING "#,W";Field
1840   GOSUB Update_field
1850   RETURN
1860 Cycle:~
1870   OFF KEY
1880   OFF KNOB
1890   GOSUB Update_field
1900   ON KEY 3 LABEL "Field",2 CALL Weird_event
1910   ON KEY 5 LABEL "Hold",4 GOTO Stop
1920   ON CYCLE 5,3 GOSUB Step
1930   WHILE True
1940   END WHILE
1950 Stop:~
1960   ON KEY 5 LABEL "Cycle",2 GOSUB Cycle
1970   ON KEY 7 LABEL "Scan",2 GOSUB Scan
1980   ON KEY 2 LABEL "Set",2 GOSUB Set
1990   ON KEY 6 LABEL "DOWN",2 GOSUB Dec
2000   ON KEY 1 LABEL "UP",2 GOSUB Inc
2010   ON KEY 0 LABEL "Step",2 GOSUB Step
2020   ON KEY 9 LABEL "Escape",2 GOTO Done
2030   ON KNOB .1,3 GOSUB Turn
2040   OFF CYCLE
2050   RETURN
2060 Update_field:~
2070   F$=VAL$(Field)
2080   ON KEY 8 LABEL F$,2 CALL Weird_event
2090   RETURN
2100 Scan:~
2110   OFF KNOB
2120   CALL Magnet_scan(X_gdu_max,Y_gdu_max,Numpeak,Interference,Mass(*),Pot(*))
2130   OUTPUT 12 USING "#,W";Pot(I)
2140   DISP USING "3D,"" = "",4Z";Mass(I);Pot(I)
2150   ON KNOB .1,3 GOSUB Turn
2160   RETURN
2170 Done:~
2180   OFF KNOB
2190   OUTPUT 12 USING "#,W";Pot(1)
2200   IF Interference THEN Pot(5)=(Pot(4)+Pot(3))/2
2210 SUBEND
2220 SUB Final_print(S$,Rcorr(*),INTEGER Ncorr,Blockmax)
2230   INTEGER N,R
2240   PRINT USING "@~"
2250   PRINT
2260   PRINT "U.S.G.S. B.I.G. #1 6""""
2270   PRINT DATE$(TIMEDATE)
2280   PRINT TIME$(TIMEDATE)
2290   PRINT TAB(20);S$
2300   PRINT
2310   FOR N=1 TO Blockmax
2320     FOR R=1 TO Ncorr
2330       PRINT USING "2X,3D.7D,#";Rcorr(R,N)
2340     NEXT R
2350     PRINT
2360     PRINT
2370   NEXT N
2380 SUBEND

```

```

2390 SUB Grandstats(R(*),Mean(*),Sigma(*),Sigma_int,INTEGER Nratio,Nblock)
2400 OPTION BASE 1
2410 IF Nblock=1 THEN SUBEXIT
2420 INTEGER N,K
2430 ALLOCATE Sumdev(Nratio),Weight(Nblock)
2440 REAL Sum_weight
2450 Sum_weight=0
2460 FOR N=1 TO Nblock
2470   Sum_weight=Sum_weight+R(7,N)
2480 NEXT N
2490 Sigma_int=1/SQR(Sum_weight)
2500 FOR N=1 TO Nblock
2510   Weight(N)=R(7,N)/Sum_weight
2520 NEXT N
2530 FOR K=1 TO Nratio
2540   Mean(K)=0
2550   FOR N=1 TO Nblock
2560     Mean(K)=Mean(K)+R(K,N)*Weight(N)
2570   NEXT N
2580   Sumdev(K)=0
2590   FOR N=1 TO Nblock
2600     Sumdev(K)=Sumdev(K)+(R(K,N)-Mean(K))^2
2610   NEXT N
2620   Sigma(K)=SQR(Sumdev(K)/(Nblock*(Nblock-1)))
2630 NEXT K
2640 SUBEND
2650 SUB Identify(Elem$,Sample$,X_gdu_max,Y_gdu_max)
2660 DISP CHR$(128);
2670 GINIT
2680 GRAPHICS ON
2690 CSIZE 50
2700 MOVE X_gdu_max/2,Y_gdu_max/2
2710 LORG 5
2720 LABEL Elem$
2730 LINPUT "Sample name (<=35 characters)?",Sample$
2740 SUBEND
2750 SUB Linear_regress(Times(*),Values(*),Cmean,Std_d,INTEGER Num,Rej)
2760 INTEGER K,J,N
2770 REAL Slope,Intcpt,Sigmult,Sumx,Sumx2,Sumy,Sumxy,Sumd,D
2780 Rej=0
2790 GOSUB Mean_sd_lr
2800 J=0
2810 Sigmult=1.16+Num*.0988-Num^2*.00214
2820 FOR K=1 TO Num
2830   D=Values(K)-Intcpt-Slope*Times(K)
2840   IF ABS(D)>Sigmult*Std_d THEN
2850     Rej=Rej+1
2860   ELSE
2870     J=J+1
2880     Values(J)=Values(K)
2890     Times(J)=Times(K)
2900   END IF
2910 NEXT K
2920 IF Rej=0 OR Num-Rej<2 THEN SUBEXIT
2930 GOSUB Mean_sd_lr
2940 SUBEXIT
2950 Mean_sd_lr:
2960 Sumx=0
2970 Sumx2=0
2980 Sumy=0
2990 Sumxy=0
3000 N=Num-Rej
3010 FOR K=1 TO N
3020   Sumx=Sumx+Times(K)
3030   Sumx2=Sumx2+Times(K)^2
3040   Sumy=Sumy+Values(K)

```

```

3050     Sumxy=Sumxy+Values(K)*Times(K)
3060 NEXT K
3070 Cmean=Sumy/N
3080 Intcpt=(Sumx2*Sumy-Sumx*Sumxy)/(N*Sumx2-Sumx^2)
3090 Slope=(N*Sumxy-Sumx*Sumy)/(N*Sumx2-Sumx^2)
3100 Sumd=0
3110 FOR K=1 TO N
3120     Sumd=Sumd+(Values(K)-Intcpt-Slope*Times(K))^2
3130 NEXT K
3140 Std_d=SQR(Sumd/(N-2))
3150 RETURN
3160 SUBEND
3170 SUB Magnet_scan(X_gdu_max,Y_gdu_max,INTEGER Numpeak,Interference,Mass(*),Pot(*))
3180     INTEGER Offset,Begin,End,Speed,True,Field,Scanup,First,Last,Tem,Logscale
3190     INTEGER Start
3200     REAL Dvm,Dplot,Del,Upper,Lower,Temp,L
3210     DIM B$(14),E$(14),F$(4),L$(6),U$(10),L1$(10),S$(10)
3220     OFF KEY
3230     ON KEY 9 LABEL "Escape",3 GOTO Done_it
3240     True=1
3250     Del=.00368*Mass(1)+.153
3260     Offset=PROUND(ABS(Pot(1)-Pot(2))/ABS(Mass(1)-Mass(2))*Del+20,0)
3270     Begin=MIN(Pot(*))-Offset
3280     End=MAX(Pot(*))+Offset
3290     GOSUB Reduce
3300     Speed=4
3310     Scanup=True
3320     Logscale=True
3330     Upper=5
3340     Lower=0
3350     GOSUB Scale
3360     GOSUB Scan_dir
3370     ON KEY 0 LABEL "Start Scan",3 GOSUB New
3380     OUTPUT 716;"ITIME USER=.5"
3390     OUTPUT 716;"HISTORY CLEAR"
3400     WHILE True
3410         ENTER 716;Dvm
3420         DISP USING "5D.3D,X,""mV""";Dvm*1000
3430     END WHILE
3440 New: !
3450     OFF KEY
3460     First=End+1
3470     IF Scanup THEN First=Begin-1
3480     Last=Begin
3490     IF Scanup THEN Last=End
3500     OUTPUT 12 USING "#,W";First
3510     GINIT
3520     CALL Clear_crt
3530     GRAPHICS OFF
3540     VIEWPORT .12*X_gdu_max,.96*X_gdu_max,.3*Y_gdu_max,.95*Y_gdu_max
3550     FRAME
3560     IF Logscale THEN
3570         WINDOW Begin,End,-1,4
3580         AXES 20,0,Begin,-1
3590         LINE TYPE 4
3600         FOR L=0 TO 3
3610             MOVE Begin,L
3620             PLOT End,L,-1
3630         NEXT L
3640         LINE TYPE 1
3650         CLIP OFF
3660         CSIZE 4
3670         LOG 6
3680         MOVE Begin,-1.05
3690         LABEL USING "4Z";Begin
3700         MOVE End,-1.05

```

```

3710 LABEL USING "4Z";End
3720 LORG 8
3730 L=-1
3740 MOVE Begin,L
3750 LABEL USING "1Z.1D,X,""mV""";10^L
3760 FOR L=0 TO 2
3770 MOVE Begin,L
3780 LABEL USING "3D,X,""mV""";10^L
3790 NEXT L
3800 FOR L=3 TO 4
3810 MOVE Begin,L
3820 LABEL USING "2D,X,""V""";10^L/1000
3830 NEXT L
3840 LORG 4
3850 FOR L=1 TO Numpeak+Interference
3860 MOVE Pot(L),4.05
3870 LABEL Mass(L)
3880 NEXT L
3890 ELSE
3900 WINDOW Begin,End,Lower,Upper
3910 AXES 20,0,Begin,Lower
3920 CLIP OFF
3930 CSIZE 4
3940 LORG 6
3950 MOVE Begin,Lower
3960 LABEL USING "4Z";Begin
3970 MOVE End,Lower
3980 LABEL USING "4Z";End
3990 LORG 8
4000 FOR L=Lower TO Upper STEP Upper-Lower
4010 MOVE Begin,L
4020 LABEL USING "2D.2D,X,""V""";L
4030 NEXT L
4040 LORG 4
4050 FOR L=1 TO Numpeak
4060 MOVE Pot(L),Upper+Upper*.002
4070 LABEL Mass(L)
4080 NEXT L
4090 END IF
4100 CLIP ON
4110 GRAPHICS ON
4120 ON KEY 3 LABEL "Field",3 CALL Weird_event
4130 ON KEY 9 LABEL "Escape",4 GOTO Done_it
4140 DISP "Waiting for magnet to stabilize..."
4150 WAIT 4
4160 OUTPUT 716;"HISTORY CLEAR"
4170 ENTER 716;Dvm
4180 DISP USING "5D.3D,X,""mV""";Dvm*1000
4190 Dplot=Dvm*1000-5
4200 IF Dplot<=0 THEN Dplot=.01
4210 IF Logscale THEN
4220 MOVE First,LGT(Dplot)
4230 ELSE
4240 MOVE First,Dplot/1000
4250 END IF
4260 Start=First+(2*Scanup-1)
4270 OUTPUT 716;"HISTORY CLEAR"
4280 ON KEY 1 LABEL "Slower",4 GOTO Slower
4290 ON KEY 6 LABEL "Faster",4 GOTO Faster
4300 Do_it:
4310 FOR Field=Start TO Last STEP Speed
4320 OUTPUT 12 USING "#,W";Field
4330 F$=VAL$(Field)
4340 ON KEY 8 LABEL F$,3 CALL Weird_event
4350 ENTER 716;Dvm
4360 DISP USING "5D.3D,X,""mV""";Dvm*1000

```

```

4370     Dplot=Dvm*1000-5
4380     IF Dplot<=0 THEN Dplot=.01
4390     IF Logscale THEN
4400         PLOT Field,LGT(Dplot),-1
4410     ELSE
4420         PLOT Field,Dplot/1000,-1
4430     END IF
4440 NEXT Field
4450 OFF KEY
4460 ON KEY 0 LABEL "Start Scan",3 GOSUB New
4470 ON KEY 2 LABEL B$,3 GOSUB Newbegin
4480 ON KEY 7 LABEL E$,3 GOSUB Newend
4490 ON KEY 5 LABEL L$,3 GOSUB Newscale
4500 IF NOT Logscale THEN
4510     ON KEY 6 LABEL U$,3 GOSUB Newmax
4520     ON KEY 1 LABEL Ll$,3 GOSUB Newmin
4530 END IF
4540 ON KEY 4 LABEL "Dump Graphics",3 CALL Dump
4550 ON KEY 9 LABEL "Escape",3 GOTO Done_it
4560 ON KEY 3 LABEL S$,3 GOSUB New_scan_dir
4570 RETURN
4580 Slower:!
4590     Speed=Speed/2
4600     IF ABS(Speed)<1 THEN Speed=1
4610     IF Scanup AND Speed<0 THEN Speed=-Speed
4620     IF NOT Scanup AND Speed>0 THEN Speed=-Speed
4630     Start=Field
4640     OUTPUT 716;"HISTORY CLEAR"
4650     GOTO Do_it
4660 Faster:!
4670     Speed=Speed*2
4680     IF ABS(Speed)>32 THEN Speed=32
4690     IF Scanup AND Speed<0 THEN Speed=-Speed
4700     IF NOT Scanup AND Speed>0 THEN Speed=-Speed
4710     Start=Field
4720     OUTPUT 716;"HISTORY CLEAR"
4730     GOTO Do_it
4740 Reduce:!
4750     IF Begin<0 THEN Begin=0
4760     IF End>4090 THEN End=4090
4770     Begin=PROUND(Begin,1)
4780     B$="LO field="&VAL$(Begin)
4790     End=PROUND(End,1)
4800     E$="HI field="&VAL$(End)
4810     ON KEY 2 LABEL B$,3 GOSUB Newbegin
4820     ON KEY 7 LABEL E$,3 GOSUB Newend
4830     RETURN
4840 Newbegin:!
4850     INPUT "New lower field value",Tem
4860     IF Tem<End THEN Begin=Tem
4870     GOSUB Reduce
4880     RETURN
4890 Newend:!
4900     INPUT "New upper field value",Tem
4910     IF Tem>Begin THEN End=Tem
4920     GOSUB Reduce
4930     RETURN
4940 Scale:!
4950     L$="LINEAR"
4960     IF Logscale THEN L$="LOG"
4970     ON KEY 5 LABEL L$,3 GOSUB Newscale
4980     IF NOT Logscale THEN
4990         U$="Max="&VAL$(Upper)&"V"
5000         Ll$="Min="&VAL$(Lower)&"V"
5010         ON KEY 6 LABEL U$,3 GOSUB Newmax
5020         ON KEY 1 LABEL Ll$,3 GOSUB Newmin

```

```

5030 ELSE
5040 OFF KEY 6
5050 OFF KEY 1
5060 END IF
5070 RETURN
5080 Newscale: !
5090 Logscale= NOT Logscale
5100 GOSUB Scale
5110 RETURN
5120 Newmin: !
5130 INPUT "New lower limit for plot (in volts)",Temp
5140 IF Temp<Upper THEN Lower=Temp
5150 GOSUB Scale
5160 RETURN
5170 Newmax: !
5180 INPUT "New upper limit for plot (in volts)",Temp
5190 IF Temp>Lower THEN Upper=Temp
5200 GOSUB Scale
5210 RETURN
5220 New_scan_dir: !
5230 Scanup= NOT Scanup
5240 Speed=-Speed
5250 GOSUB Scan_dir
5260 RETURN
5270 Scan_dir: !
5280 S$="SCAN DOWN"
5290 IF Scanup THEN S$="SCAN UP"
5300 ON KEY 3 LABEL S$,3 GOSUB New_scan_dir
5310 RETURN
5320 Done_it: !
5330 OUTPUT 716;"ITIME USER=1"
5340 SUBEND
5350 SUB Measure_inter(Block$,Baseline(*),Ipeak(*),Itime(*),INTEGER
Delay(*),Mass(*),Intn(*),Pot(*))
5360 OPTION BASE 1
5370 DIM K$[5],M$[3],F$[4]
5380 INTEGER T,P,Rej
5390 REAL Electrometer(60),Time(60),Sumtime,Cmean
5400 OFF KEY
5410 ON KEY 5 LABEL Block$ CALL Weird_event
5420 ON KEY 0 LABEL "Block" CALL Weird_event
5430 ON KEY 2 LABEL "Rubidium" CALL Weird_event
5440 ON KEY 3 LABEL "Field" CALL Weird_event
5450 P=5
5460 K$=VAL$(Mass(P))&"RB"
5470 M$=VAL$(Mass(P))
5480 F$=VAL$(Pot(P))
5490 OUTPUT 12 USING "#,W";Pot(P)
5500 ON KEY 7 LABEL M$ CALL Weird_event
5510 ON KEY 8 LABEL F$ CALL Weird_event
5520 DISP "Delay in progress, ";Rej;" readings were rejected last integration."
5530 WAIT Delay(P)
5540 Sumtime=0
5550 OUTPUT 716;"HISTORY CLEAR"
5560 FOR T=1 TO Intn(P)
5570 ENTER 716;Electrometer(T)
5580 Time(T)=TIMEDATE
5590 DISP USING "5A,""= """,5D.3D,"" mV at """,8A";K$;(Electrometer(T)-
Baseline(P,2))*1000;TIME$(Time(T))
5600 Sumtime=Sumtime+Time(T)
5610 NEXT T
5620 OUTPUT 12 USING "#,W";Pot(1)
5630 CALL Edit_routine(Electrometer(*),Cmean,(Intn(P)),Rej)
5640 Ipeak(2)=Cmean
5650 Itime(2)=Sumtime/Intn(P)
5660 SUBEND

```

```

5670 SUB Measure_peaks(Element$,Elem$,Block$,Baseline(*),Peak(*),Timep(*),INTEGER
Scan,Numpeak,Mass(*),Pot(*),Delay(*),Intn(*),Timbegin,Block,Params)
5680 OPTION BASE 1
5690 INTEGER S,P,Rej,T,True,False
5700 REAL Sumtime,Electrometer(60),Time(60),Cmean
5710 DIM K$(5),Scan$(8),M$(3),F$(4)
5720 OFF KEY
5730 False=0
5740 True= NOT False
5750 Rej=0
5760 DISP CHR$(129);
5770 ON KEY 5 LABEL Block$ CALL Weird_event
5780 ON KEY 1 LABEL "Scan" CALL Weird_event
5790 ON KEY 0 LABEL "Block" CALL Weird_event
5800 ON KEY 9 LABEL "ABORT Block" GOTO Block_abort
5810 ON KEY 2 LABEL Element$ CALL Weird_event
5820 ON KEY 3 LABEL "Field" CALL Weird_event
5830 FOR S=1 TO Scan
5840   FOR P=1 TO Numpeak
5850     K$=VAL$(Mass(P))&Elem$
5860     M$=VAL$(Mass(P))
5870     Scan$=VAL$(S)&" of "&VAL$(Scan)
5880     F$=VAL$(Pot(P))
5890     OUTPUT 12 USING "#,W";Pot(P)
5900     ON KEY 6 LABEL Scan$ CALL Weird_event
5910     ON KEY 7 LABEL M$ CALL Weird_event
5920     ON KEY 8 LABEL F$ CALL Weird_event
5930     DISP "Delay in progress, ";Rej;" readings were rejected last integration."
5940     WAIT Delay(P)
5950     Sumtime=0
5960     OUTPUT 716;"HISTORY CLEAR"
5970     FOR T=1 TO Intn(P)
5980       ENTER 716;Electrometer(T)
5990       Time(T)=TIMEDATE
6000       DISP USING "5A,""= """,5D.3D,"" mV at """,8A";K$(Electrometer(T)-
Baseline(P,1))*1000;TIME$(Time(T))
6010       Sumtime=Sumtime+Time(T)
6020     NEXT T
6030     CALL Edit_routine(Electrometer(*),Cmean,(Intn(P)),Rej)
6040     Peak(P,S)=Cmean
6050     Timep(P,S)=Sumtime/Intn(P)
6060     IF S=1 THEN
6070       MOVE ((Timep(P,S) MOD 86400)/60-Timbegin)-(Intn(P)/2)/60,(Peak(P,S)-
Baseline(P,1))
6080       GOTO First_scan
6090     END IF
6100     MOVE ((Timep(P,S-1) MOD 86400)/60-Timbegin)+(Intn(P)/2)/60,(Peak(P,S-1)-
Baseline(P,1))
6110     LINE TYPE 4
6120     PLOT ((Timep(P,S) MOD 86400)/60-Timbegin)-(Intn(P)/2)/60,(Peak(P,S)-
Baseline(P,1)), -1
6130 First_scan:
6140     LINE TYPE 1
6150     IDRAW Intn(P)/60,0
6160     NEXT P
6170   NEXT S
6180 SUBEXIT
6190 Block_abort:
6200   Params=True
6210   Block=Block-1
6220 SUBEND
6230 SUB Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),INTEGER
Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
6240 OPTION BASE 1
6250 INTEGER P,S,Rej,T,B_o,First,Last,Offset,Field
6260 REAL Sumtime,Electrometer(120),Time(120),Zero(5,2),Timezero(5,2),Cmean

```

```

6270 REAL Del
6280 DIM M$(3),K$(5),F$(4)
6290 OFF KEY
6300 DISP CHR$(129);
6310 ON KEY 0 LABEL "Block" CALL Weird_event
6320 ON KEY 5 LABEL Block$ CALL Weird_event
6330 ON KEY 7 LABEL "Baselines" CALL Weird_event
6340 ON KEY 2 LABEL Element$ CALL Weird_event
6350 ON KEY 3 LABEL "Field" CALL Weird_event
6360 Del=.00368*Mass(1)+.153
6370 Offset=PROUND(ABS(Pot(1)-Pot(2))/ABS(Mass(1)-Mass(2))*Del,0)
6380 IF Mass(1)<=Mass(2) THEN
6390     First=2
6400     Last=1
6410 ELSE
6420     First=1
6430     Last=2
6440 END IF
6450 FOR P=Numpeak+Interference TO 1 STEP -1
6460     FOR S=First TO Last STEP Last-First
6470         B_o=(S-1)*Offset+(S-2)*Offset
6480         M$=VAL$(Mass(P))
6490         IF S=2 THEN
6500             K$="Above"
6510         ELSE
6520             K$="Below"
6530         END IF
6540         Field=Pot(P)+B_o
6550         F$=VAL$(Field)
6560         OUTPUT 12 USING "#,W";Field
6570         ON KEY 4 LABEL K$ CALL Weird_event
6580         ON KEY 9 LABEL M$ CALL Weird_event
6590         ON KEY 8 LABEL F$ CALL Weird_event
6600         DISP "Delay in progress, ";Rej;" readings were rejected last integration."
6610         WAIT Delay(1)
6620         Sumtime=0
6630         OUTPUT 716;"HISTORY CLEAR"
6640         FOR T=1 TO Intn(P)*2
6650             ENTER 716;Electrometer(T)
6660             Time(T)=TIMEDATE
6670             DISP USING """"Baseline= "",5D.3D,"" mV at
6680             """,8A";Electrometer(T)*1000;TIME$(Time(T))
6690             Sumtime=Sumtime+Time(T)
6700         NEXT T
6710         CALL Edit_routine(Electrometer(*),Cmean,Intn(P)*2,Rej)
6720         Zero(P,S)=Cmean
6730         Timezero(P,S)=Sumtime/(Intn(P)*2)
6740     NEXT S
6750 NEXT P
6760 IF NOT Interference THEN OUTPUT 12 USING "#,W";Pot(1)
6770 FOR P=1 TO Numpeak+Interference
6780     Baseline(P,2)=(Zero(P,1)+Zero(P,2))/2
6790     Timebase(P,2)=(Timezero(P,1)+Timezero(P,2))/2
6800 NEXT P
6810 SUB New_zeros(Baseline(*),Timebase(*),INTEGER Numpeak,Interference,OPTIONAL
Ipeak(*),Itime(*)
6820     INTEGER P
6830     IF Interference THEN
6840         Ipeak(1)=Ipeak(2)
6850         Itime(1)=Itime(2)
6860     END IF
6870     FOR P=1 TO Numpeak+Interference
6880         Baseline(P,1)=Baseline(P,2)
6890         Timebase(P,1)=Timebase(P,2)
6900     NEXT P

```

```

6910 SUBEND
6920 SUB Parameters (Elem$, INTEGER
Check, Numpeak, Interference, Blockmax, Scan, Factor, Mass (*), Intn (*), Delay (*), OPTIONAL Disc)
6930   INTEGER N, True, False
6940   DIM Ans1$(3), Ans2$(3), Ans3$(3)
6950   CALL Clear_crt
6960   OFF KEY
6970   False=0
6980   True= NOT False
6990   DISP CHR$(128);
7000   OUTPUT 1; "                                CURRENT RUN PARAMETERS"
7010   OUTPUT 1
7020   OUTPUT 1; "Total number of blocks="; Blockmax
7030   OUTPUT 1
7040   OUTPUT 1; "Number of scans per block="; Scan
7050   OUTPUT 1
7060   OUTPUT 1; "Integration times (seconds):"
7070   OUTPUT 1; " ";
7080   FOR N=1 TO Numpeak+Interference
7090     OUTPUT 1; "   "; Mass(N); "="; Intn(N);
7100   NEXT N
7110   OUTPUT 1 USING "/"
7120   OUTPUT 1; "Delay times (seconds):"
7130   OUTPUT 1; " ";
7140   FOR N=1 TO Numpeak+Interference
7150     OUTPUT 1; "  Before"; Mass(N); "="; Delay(N);
7160   NEXT N
7170   OUTPUT 1 USING "/"
7180   IF Factor THEN OUTPUT 1; "Complete data output."
7190   IF NOT Factor THEN OUTPUT 1; "Print means only."
7200   OUTPUT 1
7210   IF Interference AND Elem$="SR" AND SIZE(Mass,1)=5 THEN OUTPUT 1; "Correct for
87Rb."
7220   IF NOT Interference AND Elem$="SR" AND SIZE(Mass,1)=5 THEN OUTPUT 1; "No Rb
correction."
7230   IF NPAR>10 THEN OUTPUT 1; "Mass discrimination (per amu) =" ; Disc
7240   INPUT "Do you want to change any of these parameters?", Ans1$
7250   IF UPC$(Ans1$)="N" OR UPC$(Ans1$)="NO" THEN
7260     Check=True
7270     SUBEXIT
7280   END IF
7290   INPUT "Number of blocks (<=50)", Blockmax
7300   INPUT "Number of scans per block (<=25)", Scan
7310   DISP "Integration times in sec (" ;
7320   FOR N=1 TO Numpeak+Interference
7330     DISP Mass(N); ", ";
7340   NEXT N
7350   DISP CHR$(8); ")";
7360   INPUT Intn(*)
7370   DISP "Delay times in sec (" ;
7380   FOR N=1 TO Numpeak+Interference
7390     DISP Mass(N); ", ";
7400   NEXT N
7410   DISP CHR$(8); ")";
7420   INPUT Delay(*)
7430   INPUT "Complete data output (Yes or No)", Ans2$
7440   IF UPC$(Ans2$)="Y" OR UPC$(Ans2$)="YES" THEN
7450     Factor=True
7460   ELSE
7470     Factor=False
7480   END IF
7490   IF Elem$="SR" AND SIZE(Mass,1)=5 THEN
7500     INPUT "Correct for 87Rb interference", Ans3$
7510     IF UPC$(Ans3$)="N" OR UPC$(Ans3$)="NO" THEN
7520       Interference=False
7530     ELSE

```

```

7540     Interference=True
7550     END IF
7560     END IF
7570     IF NPAR>10 THEN
7580         INPUT "Mass discrimination (per amu)",Disc
7590     END IF
7600     Check=False
7610 SUBEND
7620 SUB Plot (Qpeak(*),X_gdu_max,Y_gdu_max,INTEGER Timfirst,Timlast)
7630     REAL Maxplot,L
7640     SELECT MAX(Qpeak(*))*1.3
7650     CASE <.1
7660         Maxplot=.1
7670     CASE <1
7680         Maxplot=1
7690     CASE ELSE
7700         Maxplot=10
7710     END SELECT
7720     GINIT
7730     CALL Clear_crt
7740     GRAPHICS OFF
7750     VIEWPORT .15*X_gdu_max,.99*X_gdu_max,.25*Y_gdu_max,.9*Y_gdu_max
7760     FRAME
7770     WINDOW Timfirst,Timlast,0,Maxplot
7780     AXES .1,Maxplot/50,Timfirst,Maxplot,10,5,4
7790     CLIP OFF
7800     DEG
7810     LDIR 90
7820     MOVE Timfirst-(Timlast-Timfirst)*.12,Maxplot/2
7830     LORG 4
7840     LABEL "INTENSITY (V)"
7850     LDIR 0
7860     MOVE Timfirst+(Timlast-Timfirst)/2,Maxplot+Maxplot*.04
7870     LABEL "Time since start of run (minutes)"
7880     CSIZE 4
7890     FOR L=Timfirst TO Timlast
7900         MOVE L,Maxplot+Maxplot*.002
7910         LABEL L
7920     NEXT L
7930     LORG 8
7940     FOR L=0 TO Maxplot STEP Maxplot/10
7950         MOVE Timfirst-.05,L
7960         LABEL L
7970     NEXT L
7980     CLIP ON
7990     GRAPHICS ON
8000 SUBEND
8010 SUB Print_means (Mean(*),Sigma(*),Sigma_int,INTEGER Nblock,Nratio)
8020     INTEGER R
8030     IF Nblock=1 THEN
8040         PRINT USING "6(/)"
8050         SUBEXIT
8060     END IF
8070     PRINT
8080     PRINT "Weighted means and external 2-sigmas for ";Nblock;" blocks:"
8090     FOR R=1 TO Nratio
8100         PRINT USING "4X, 3D.7D, #";Mean(R)
8110     NEXT R
8120     PRINT
8130     FOR R=1 TO Nratio
8140         PRINT USING "4X, 3D.7D, #";2*Sigma(R)
8150     NEXT R
8160     PRINT USING "4X, 3D.6D";2*Sigma_int
8170     PRINT USING "6(/)"
8180 SUBEND
8190 SUB Print_out1 (Traw$(*),Baseline(*),Netpeak(*),Rraw(*),INTEGER

```

```

Factor, Numpeak, Interference, Block, Scan, Mass (*), Pot (*)
8200  OPTION BASE 1
8210  INTEGER N, K
8220  DIM M$(4) [7]
8230  PRINT "Block "; Block; " finished at "; TIME$(TIMEDATE)
8240  FOR N=1 TO Numpeak+Interference
8250    PRINT TAB(32+(N-1)*10); Mass(N);
8260  NEXT N
8270  PRINT
8280  PRINT "Baselines before this block:";
8290  FOR N=1 TO Numpeak+Interference
8300    PRINT USING "X, 5D.3D, #"; Baseline(N, 1)*1000
8310  NEXT N
8320  PRINT
8330  PRINT "Baselines after this block: ";
8340  FOR N=1 TO Numpeak+Interference
8350    PRINT USING "X, 5D.3D, #"; Baseline(N, 2)*1000
8360  NEXT N
8370  PRINT
8380  PRINT "Initial peak intensities: ";
8390  FOR N=1 TO Numpeak+Interference
8400    PRINT USING "X, 5D.3D, #"; Netpeak(N, 1)*1000
8410  NEXT N
8420  PRINT
8430  PRINT "Final peak intensities: ";
8440  FOR N=1 TO Numpeak+Interference
8450    PRINT USING "X, 5D.3D, #"; Netpeak(N, Scan)*1000
8460  NEXT N
8470  PRINT
8480  PRINT "Magnet field values: ";
8490  FOR N=1 TO Numpeak+Interference
8500    PRINT USING "2X, 4Z, 4X, #"; Pot(N)
8510  NEXT N
8520  PRINT
8530  PRINT
8540  IF NOT Factor THEN SUBEXIT
8550  FOR N=2 TO Numpeak
8560    M$(N) = VAL$(Mass(N)) & "/" & VAL$(Mass(1))
8570  NEXT N
8580  FOR N=2 TO Numpeak
8590    PRINT USING "4X, 7A, 5X, ""Time"", #"; M$(N)
8600  NEXT N
8610  PRINT
8620  FOR K=1 TO Scan-1
8630    FOR N=2 TO Numpeak
8640      PRINT USING "3D.6D, 2X, 8A, #"; Rraw(N, K); Traw$(N, K)
8650    NEXT N
8660    PRINT
8670  NEXT K
8680 SUBEND
8690 SUB Print_out2(Meanraw(*), Percent_dev(*), INTEGER Numpeak, Rejected(*), Mass(*))
8700  OPTION BASE 1
8710  DIM M$(4) [7]
8720  INTEGER N
8730  FOR N=2 TO Numpeak
8740    M$(N) = VAL$(Mass(N)) & "/" & VAL$(Mass(1))
8750  NEXT N
8760  FOR N=2 TO Numpeak
8770    PRINT USING "4X, 7A, 4X, ""% Dev"", #"; M$(N)
8780  NEXT N
8790  PRINT
8800  FOR N=2 TO Numpeak
8810    PRINT USING "3D.7D, 3D.5D, #"; Meanraw(N); Percent_dev(N)
8820  NEXT N
8830  PRINT
8840  IF MAX(Rejected(*))=0 THEN

```

```

8850     PRINT "No ratios were rejected."
8860 ELSE
8870     PRINT "Rejected:";
8880     FOR N=2 TO Numpeak
8890         PRINT TAB(15+(N-2)*20);Rejected(N);
8900     NEXT N
8910     PRINT
8920 END IF
8930 SUBEND
8940 SUB
Ratios(Traw$(*),Baseline(*),Timebase(*),Timep(*),Peak(*),Netpeak(*),Rraw(*),Avetime(*),INT
EGER Numpeak,Scan,Interference,OPTIONAL Ipeak(*),Itime(*))
8950     OPTION BASE 1
8960     INTEGER S,P,Numb
8970     REAL Zero_rate(5),Base(5,25),Peak_rate(5,25),Irate,Ibase
8980     FOR P=1 TO Numpeak+Interference
8990         Zero_rate(P)=(Baseline(P,2)-Baseline(P,1))/(Timebase(P,2)-Timebase(P,1))
9000     NEXT P
9010     FOR P=1 TO Numpeak
9020         FOR S=1 TO Scan
9030             Base(P,S)=Baseline(P,1)+Zero_rate(P)*(Timep(P,S)-Timebase(P,1))
9040             Netpeak(P,S)=Peak(P,S)-Base(P,S)
9050         NEXT S
9060     NEXT P
9070     IF Interference THEN
9080         FOR S=1 TO 2
9090             Base(5,S)=Baseline(5,1)+Zero_rate(5)*(Itime(S)-Timebase(5,1))
9100         NEXT S
9110         Netpeak(5,1)=Ipeak(1)-Base(5,1)
9120         Netpeak(5,Scan)=Ipeak(2)-Base(5,2)
9130         Irate=(Netpeak(5,Scan)-Netpeak(5,1))/(Itime(2)-Itime(1))
9140         FOR S=1 TO Scan
9150             Ibase=Netpeak(5,1)+Irate*(Timep(2,S)-Itime(1))
9160             IF Ibase>0 THEN Netpeak(2,S)=Netpeak(2,S)-Ibase*.387
9170         NEXT S
9180     END IF
9190     Numb=Scan-1
9200     FOR P=1 TO Numpeak
9210         FOR S=1 TO Numb
9220             Peak_rate(P,S)=(Netpeak(P,S)-Netpeak(P,S+1))/(Timep(P,S)-Timep(P,S+1))
9230         NEXT S
9240     NEXT P
9250     FOR P=2 TO Numpeak
9260         FOR S=1 TO Numb
9270             Avetime(P,S)=(Timep(P,S)+Timep(P,S+1)+Timep(1,S)+Timep(1,S+1))/4
9280             Rraw(P,S)=(Netpeak(P,S)+Peak_rate(P,S)*(Avetime(P,S)-
Timep(P,S)))/(Netpeak(1,S)+Peak_rate(1,S)*(Avetime(P,S)-Timep(1,S)))
9290             Traw$(P,S)=TIME$(Avetime(P,S))
9300         NEXT S
9310     NEXT P
9320 SUBEND
9330 SUB Start_up(Block$,Sample$,INTEGER Block,Blockmax,Timbegin)
9340     CALL Clear_crt
9350     DISP CHR$(128);
9360     DISP TAB(20);CHR$(131);"Printer ON LINE?"
9370     WAIT 5
9380     DISP CHR$(128);CHR$(12);
9390     PRINT TAB(20);Sample$;TAB(60);DATE$(TIMEDATE)
9400     PRINT
9410     Timbegin=INT((TIMEDATE MOD 86400)/60)
9420     Block=1
9430     Block$=VAL$(Block)&" of "&VAL$(Blockmax)
9440 SUBEND
9450 SUB Statistics(Avetime(*),Rraw(*),Meanraw(*),Percent_dev(*),INTEGER
Numpeak,Scan,Rejected(*))
9460     OPTION BASE 1

```

```

9470  INTEGER P,K,Numb,Rej
9480  REAL Regress_time,Ratio(24),Tim(24),Cmean,Std_d,Std_dev
9490  Numb=Scan-1
9500  Regress_time=TIMEDATE
9510  FOR P=2 TO Numpeak
9520    FOR K=1 TO Numb
9530      Ratio(K)=Rraw(P,K)
9540      Tim(K)=Avetime(P,K)-Regress_time
9550    NEXT K
9560    CALL Linear_regress(Tim(*),Ratio(*),Cmean,Std_d,(Numb),Rej)
9570    Rejected(P)=Rej
9580    Meanraw(P)=Cmean
9590    Std_dev=Std_d/SQR(Numb-Rejected(P))
9600    Percent_dev(P)=Std_dev/Meanraw(P)*100
9610  NEXT P
9620  SUBEND
9630  SUB Summ_plot(Rcorr(*),Mean(*),Sigma(*),X_gdu_max,Y_gdu_max,INTEGER
Blockmax,Index1,Index2)
9640    OPTION BASE 1
9650    INTEGER I,L,True
9660    REAL Minplot,Maxplot,Plot_mean,Plot_sdev
9670    ALLOCATE Plot_corr(Blockmax),Plot_sigma(Blockmax)
9680    True=1
9690    FOR I=1 TO Blockmax
9700      Plot_corr(I)=Rcorr(Index1,I)
9710      Plot_sigma(I)=Rcorr(Index2,I)*Plot_corr(I)*2/100
9720    NEXT I
9730    Plot_mean=Mean(Index1)
9740    Plot_sdev=Sigma(Index1)
9750    GINIT
9760    CALL Clear crt
9770    GRAPHICS OFF
9780    VIEWPORT .01*X_gdu_max,.99*X_gdu_max,.3*Y_gdu_max,.99*Y_gdu_max
9790    FRAME
9800    Minplot=MIN(Plot_corr(*))-MAX(Plot_sigma(*))
9810    Maxplot=MAX(Plot_corr(*))+MAX(Plot_sigma(*))
9820    WINDOW 0,Blockmax+1,Minplot,Maxplot
9830    AXES 1,0,0,Minplot,1,1,4
9840    CLIP OFF
9850    CSIZE 4
9860    LORG 6
9870    FOR L=5 TO Blockmax+1 STEP 5
9880      MOVE L,Minplot
9890      LABEL L
9900    NEXT L
9910    CLIP ON
9920    FOR I=1 TO Blockmax
9930      MOVE I-.3,Plot_corr(I)-Plot_sigma(I)
9940      RECTANGLE .6,2*Plot_sigma(I),FILL,EDGE
9950    NEXT I
9960    FOR I=-1 TO 1
9970      MOVE 0,Plot_mean+I*2*Plot_sdev
9980      LINE TYPE ABS(I)*3+1
9990      IDRAW Blockmax+1,0
10000  NEXT I
10010  GRAPHICS ON
10020  ON KEY 4 LABEL "Dump Graphics" GOTO Dump
10030  ON KEY 0 LABEL "Continue" GOTO Done
10040  WHILE True
10050    END WHILE
10060  Dump: !
10070    CALL Dump
10080  Done: !
10090  SUBEND
10100  SUB Weird_event
10110    BEEP

```

10120 SUBEND

KBDM

```

10   REM KBDM                      Brian D. Marshall 7/88
20   OPTION BASE 1
30   PRINTER IS PRT
40   CONTROL 1,4;0                !Turns off DISPLAY FUNCTIONS
50   LOADSUB ALL FROM "Subfile"
60   DIM
A$[4],Block$[8],Elem$[2],Element$[10],File$[6],N$[6],Sample$[35],Traw$(2,24)[8],Fil$[1]
70   INTEGER Block,Blockmax,Check,Delay(2),Edit1,Factor,False,Halt
80   INTEGER Index1,Index2,Interference,Intn(2),Mass(2),Numpeak,Params,Pot(2)
90   INTEGER Prog,Rejected(2),Scan,Timbegin,Timfirst,Timlast,True
100  REAL Avetime(2,24),Baseline(2,2),Disc,Mean(2)
110  REAL Meanraw(2),Netpeak(2,24),Peak(2,24),Percent_dev(2)
120  REAL Qpeak(2),Rcorr(7,50),Rraw(2,24),Rspike,Rstore(4),Scantime
130  REAL Sigma(2),Sigma_int,Tcorr(2),Timebase(2,2),Timep(2,24)
140  REAL X_gdu_max,Y_gdu_max
150  CALL Clear crt
160  Y_gdu_max=100*MAX(1,1/RATIO)
170  X_gdu_max=100*MAX(1,RATIO)
180  False=0
190  True= NOT False
200  Element$="Potassium"
210  Elem$="K"
220  CALL Identify(Elem$,Sample$,X_gdu_max,Y_gdu_max)
230  INPUT "Triple filament as perchlorate (T) or single filament as chloride (S)?",Fil$
240  ON KEY 0 LABEL "Sample" GOTO Unspiked
250  ON KEY 2 LABEL "Spiked sample" GOTO Spiked
260  ON KEY 4 LABEL "Spike" GOTO Spike
270  WHILE True
280  END WHILE
290  Unspiked:!
300  Prog=1
310  GOTO Initialize
320  Spiked:!
330  Prog=3
340  GOTO Initialize
350  Spike:!
360  Prog=2
370  Initialize:!
380  OFF KEY
390  GRAPHICS OFF
400  IF Prog=3 THEN GOSUB Spikeratios
410  GOSUB Defaults
420  ON KEY 2 LABEL "START" GOTO Main
430  ON KEY 9 LABEL "Escape" GOTO Initialize
440  ON KEY 5 LABEL "ABORT RUN!" GOTO Escape
450  WHILE True
460  END WHILE
470  Massdisc:!
480  Rcorr(1,Block)=1
490  Rcorr(2,Block)=Meanraw(2)*(1+2*Disc)
500  Rcorr(3,Block)=Percent_dev(2)
510  Rcorr(4,Block)=Block
520  Rcorr(7,Block)=1/Rcorr(3,Block)^2
530  RETURN
540  Spikesub:!
550  Tcorr(2)=Meanraw(2)*(1+2*Disc)
560  Rcorr(1,Block)=(1-Tcorr(2)*Rspike)/(Tcorr(2)-.072168)
570  Rcorr(2,Block)=(Tcorr(2)*Rspike-1)/Rcorr(1,Block)+Tcorr(2)
580  Rcorr(3,Block)=Percent_dev(2)
590  Rcorr(4,Block)=Block
600  Rcorr(7,Block)=1/Rcorr(3,Block)^2
610  RETURN
620  Printout3: !

```

!39S/41T
!41/39

```

630 Format3: IMAGE 2(4X,3D.7D)
640 IF Prog=3 THEN
650   PRINT
660   PRINT "Spike ratio: ";DROUND(Rspike,6)
670 END IF
680 PRINT
690 PRINT "Mass discrimination corrected (=";Disc;"per a.m.u.)."
700 PRINT
710 PRINT USING "2(9X,7A)";"39S/41T";" 41/39 "
720 PRINT USING Format3;Rcorr(1,Block);Rcorr(2,Block)
730 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Block),2)
740 RETURN
750 Defaults: !
760 Blockmax=10
770 Intn(1)=5
780 Intn(2)=5
790 Delay(1)=5
800 Delay(2)=5
810 Pot(1)=1250
820 Pot(2)=2170
830 Mass(1)=39
840 Mass(2)=41
850 Scan=11
860 Factor=False
870 Interference=False
880 Numpeak=2
890 IF UPC$(Fil$)="T" THEN
900   Disc=.003
910 ELSE
920   Disc=.012
930 END IF
940 Showdefaults: !
950 CALL Field_values(X_gdu_max,Y_gdu_max,Numpeak,Interference,Mass(*),Pot(*))
960 CALL
Parameters(Elem$,Check,Numpeak,Interference,Blockmax,Scan,Factor,Mass(*),Intn(*),Delay(*),
Disc)
970 IF Check THEN RETURN
980 IF NOT Check THEN GOTO Showdefaults
990 Spikeratios: !
1000 Rspike=.02015      !39/41
1010 RETURN
1020 REM Key_handlers
1030 Pause: !
1040 OFF KEY
1050 Halt=True
1060 RETURN
1070 Last_block: !
1080 OFF KEY
1090 Blockmax=Block
1100 RETURN
1110 Escape: !
1120 OFF KEY
1130 CALL Clear_crt
1140 DISP CHR$(131);"PROGRAM ABORTED"
1150 WAIT 5
1160 GOTO Lastline
1170 New_params: !
1180 OFF KEY
1190 Params=True
1200 RETURN
1210 Summary: !
1220 Index1=(Prog<>3)+1
1230 Index2=3
1240 CALL Final_print(Sample$,Rcorr(*),4,Blockmax)
1250 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Blockmax),2)
1260 CALL Summ_plot(Rcorr(*),Mean(*),Sigma(*),X_gdu_max,Y_gdu_max,Blockmax,Index1,Index2)

```

```

1270 DISP CHR$(128)
1280 LOOP
1290   Edit1=0
1300   INPUT "Number of blocks to be edited",Edit1
1310   EXIT IF Edit1=0
1320   CALL Edit(Rcorr(*),Mean(*),Sigma(*),Sigma_int,(Edit1),2,Blockmax)
1330 END LOOP
1340 RETURN
1350 Main:
1360 OFF KEY
1370 CALL Start_up(Block$,Sample$,Block,Blockmax,Timbegin)
1380 CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
1390 CALL New_zeros(Baseline(*),Timebase(*),Numpeak,Interference)
1400 WHILE Block<=Blockmax
1410   OFF KEY
1420   Halt=False
1430   Params=False
1440   CALL Center_peaks(Qpeak(*),Numpeak,Interference,Mass(*),Pot(*))
1450   Timfirst=INT((TIMEDATE MOD 86400)/60)-Timbegin
1460   Scantime=(SUM(Delay)+SUM(Intn))/60
1470   Timlast=Timfirst+PROUND(Scantime*(Scan+4),0)
1480   CALL Plot(Qpeak(*),X_gdu_max,Y_gdu_max,Timfirst,Timlast)
1490   Block$=VAL$(Block)&" of "&VAL$(Blockmax)
1500   CALL
Measure_peaks(Element$,Elem$,Block$,Baseline(*),Peak(*),Timep(*),Scan,Numpeak,Mass(*),Pot(*),Delay(*),Intn(*),Timbegin,Block,Params)
1510   IF Params THEN GOTO Hold
1520   CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
1530   ON KEY 9 LABEL "ABORT RUN!" GOTO Escape
1540   ON KEY 5 LABEL "FINISH" GOSUB Last_block
1550   ON KEY 2 LABEL "PAUSE" GOSUB Pause
1560   ON KEY 7 LABEL "Change Params" GOSUB New_params
1570   BEEP 880,2
1580   DISP CHR$(128);"Calculating results"
1590   CALL
Ratios(Traw$(*),Baseline(*),Timebase(*),Timep(*),Peak(*),Netpeak(*),Rraw(*),Avetime(*),Numpeak,Scan,Interference)
1600   CALL
Print_out1(Traw$(*),Baseline(*),Netpeak(*),Rraw(*),Factor,Numpeak,Interference,Block,Scan,Mass(*),Pot(*))
1610   CALL
Statistics(Avetime(*),Rraw(*),Meanraw(*),Percent_dev(*),Numpeak,Scan,Rejected(*))
1620   CALL Print_out2(Meanraw(*),Percent_dev(*),Numpeak,Rejected(*),Mass(*))
1630   IF Prog=3 THEN
1640     GOSUB Spikesub
1650   ELSE
1660     GOSUB Massdisc
1670   END IF
1680   CALL Grandstats(Rcorr(*),Mean(*),Sigma(*),Sigma_int,2,(Block))
1690   GOSUB Printout3
1700 Hold:
1710   IF Halt THEN
1720     DISP CHR$(128);"Press CONTINUE to resume data acquisition."
1730     PAUSE
1740   END IF
1750   IF Params THEN
1760     CALL Clear_crt
1770     GOSUB Showdefaults
1780     DISP CHR$(128);"Press CONTINUE to resume data acquisition."
1790     PAUSE
1800   END IF
1810   Block=Block+1

```

```

1820   Block$=VAL$(Block)&" of "&VAL$(Blockmax)
1830   IF Params AND Block<=Blockmax THEN CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*
),Intn(*),Pot(*))
1840   CALL New_zeros(Baseline(*),Timebase(*),Numpeak,Interference)
1850 END WHILE
1860 OFF KEY
1870 GOSUB Summary
1880 Lastline:~
1890 CALL Clear crt
1900 PRINT USING "@~"
1910 LOAD "CONTROL"
1920 END

```

CABDM

```

10   REM CABDM                      Brian D. Marshall 7/88
20   OPTION BASE 1
30   PRINTER IS PRT
40   CONTROL 1,4;0                  !Turns off DISPLAY FUNCTIONS.
50   LOADSUB ALL FROM "Subfile"
60   DIM
AS[4],Block$[8],Elem$[2],Element$[10],File$[6],N$[6],Sample$[35],Traw$(4,24)[8],Today$[11]
,Now$[8],Mon$[3]
70   INTEGER B,Block,Blockmax,Check,Delay(4),Edit1,Factor,False,Halt,I
80   INTEGER Index1,Index2,Interference,Intn(4),Iter,Mass(4),Numpeak,Params
90   INTEGER Pot(4),Prog,Rejected(4),Scan,Sp,Timbegin,Timfirst,Timlast,True
100  REAL Alpha,Amu(4),Avetime(4,24),Baseline(4,2),Beta,Disc,Mean(4)
110  REAL Meancorr(4),Meanraw(4),Netpeak(4,24),P_exp,Peak(4,24),Percent_dev(4)
120  REAL Qpeak(4),Rcorr(7,50),Rraw(4,24),Rspike(3),Rstore(6),Scantime
130  REAL Sigma(4),Sigma_int,Sum_p,Tcorr(3),Timebase(4,2),Timep(4,24)
140  REAL X_gdu_max,Y_gdu_max
150  CALL Clear crt
160  Y_gdu_max=100*MAX(1,1/RATIO)
170  X_gdu_max=100*MAX(1,RATIO)
180  False=0
190  True= NOT False
200  Elem$="CA"
210  Element$="Calcium"
220  OFF KEY
230  CALL Identify(Elem$,Sample$,X_gdu_max,Y_gdu_max)
250  ON KEY 0 LABEL "Sample" GOTO Unspiked
260  ON KEY 2 LABEL "Spiked sample" GOTO Spiked
270  ON KEY 4 LABEL "Spike" GOTO Spike
280  WHILE True
290  END WHILE
300  Unspiked:~
310  Prog=1
320  GOTO Initialize
330  Spiked:~
340  Prog=3
350  GOTO Initialize
360  Spike:~
370  Prog=2
380  Initialize:~
390  OFF KEY
400  GRAPHICS OFF
410  IF Prog=3 THEN GOSUB Spikeratios
420  GOSUB Defaults
430  ON KEY 2 LABEL "START" GOTO Main
440  ON KEY 9 LABEL "Escape" GOTO Initialize
450  ON KEY 5 LABEL "ABORT RUN!" GOTO Escape
460  WHILE True
470  END WHILE
480  Massdisc:~
490  Sum_p=LOG(.31221/(Meanraw(2)/Meanraw(4)))/LOG(Amu(2)/Amu(4))

```

```

500 IF Prog=2 THEN Sum_p=0 !No correction for spike run
510 Rcorr(1,Block)=1 !42S/43T
520 Rcorr(2,Block)=1/(Meanraw(2))*(Amu(1)/Amu(2))^Sum_p !40/42
530 Rcorr(3,Block)=1/(Meanraw(4))*(Amu(1)/Amu(4))^Sum_p !40/44
540 Rcorr(4,Block)=(Meanraw(3)/Meanraw(2))*(Amu(3)/Amu(2))^Sum_p !43/42
550 Rcorr(5,Block)=SQRT(2*Percent_dev(2)^2+Percent_dev(4)^2)
560 Rcorr(6,Block)=Block
570 Rcorr(7,Block)=1/Rcorr(5,Block)^2
580 RETURN
590 Spikesub:!
600 FOR I=2 TO 4
610 Meancorr(I)=Meanraw(I)
620 NEXT I
630 Sum_p=0
640 Iter=0
650 REPEAT
660 Alpha=(Meancorr(3)/Meancorr(2)-.207745)/(1-Meancorr(3)/Meancorr(2)*Rspike(1))
670 Beta=1/(Meancorr(4)/Meancorr(2)+Alpha*Rspike(1)*(Meancorr(4)/Meancorr(2)-
Rspike(3)/Rspike(2)))
680 P_exp=LOG(.31221/Beta)/LOG(Amu(2)/Amu(4))
690 Tcorr(1)=1/Alpha
700 Tcorr(2)=1/(Meancorr(2))*(Amu(1)/Amu(2))^P_exp
710 Tcorr(3)=1/(Meancorr(4))*(Amu(1)/Amu(4))^P_exp
720 Sum_p=Sum_p+P_exp
730 Iter=Iter+1
740 Meancorr(2)=1/Tcorr(2)
750 Meancorr(3)=Meancorr(3)*(Amu(3)/Amu(1))^P_exp
760 Meancorr(4)=1/Tcorr(3)
770 UNTIL ABS(P_exp)<1.E-4
780 DISP "Number of iterations was ";Iter
790 Rcorr(1,Block)=Tcorr(1)
800 Rcorr(2,Block)=Tcorr(2)+Alpha*Rspike(1)*(Tcorr(2)-1/Rspike(2))
810 Rcorr(3,Block)=Rcorr(2,Block)*.31221
820 Rcorr(4,Block)=.207745
830 Rcorr(5,Block)=SQRT(2*Percent_dev(2)^2+Percent_dev(4)^2)
840 Rcorr(6,Block)=Block
850 Rcorr(7,Block)=1/Rcorr(5,Block)^2
860 RETURN
870 Printout3: !
880 Disc=((Amu(2)/Amu(4))^Sum_p-1)/2
890 Format3: IMAGE 5(4X,3D.7D)
900 IF Prog=3 THEN
910 PRINT
920 PRINT "Spike ratios:
";DROUND(Rspike(1),6);",";DROUND(Rspike(2),6);",";DROUND(Rspike(3),6)
930 PRINT "Number of iterations was ";Iter
940 END IF
950 PRINT
960 PRINT "Mass discrimination corrected (p=";DROUND(Sum_p,6);",
alpha=";DROUND(Disc,6);")."
970 PRINT
980 PRINT USING "5(8X,7A)";"42S/43T";" 40/42 ";" 40/44 ";" 43/42 ";"% Error"
990 PRINT USING
Format3;Rcorr(1,Block);Rcorr(2,Block);Rcorr(3,Block);Rcorr(4,Block);Rcorr(5,Block)
1000 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Block),4)
1010 RETURN
1020 Defaults:!
1030 Blockmax=30
1040 Intn(1)=5
1050 Intn(2)=9
1060 Intn(3)=7
1070 Intn(4)=7
1080 Delay(1)=5
1090 Delay(2)=6
1100 Delay(3)=4
1110 Delay(4)=4

```

```

1120 Pot(1)=1700
1130 Pot(2)=2600
1140 Pot(3)=3040
1150 Pot(4)=3475
1160 Mass(1)=40
1170 Mass(2)=42
1180 Mass(3)=43
1190 Mass(4)=44
1200 Amu(1)=39.962591
1210 Amu(2)=41.958622
1220 Amu(3)=42.958770
1230 Amu(4)=43.955485
1240 Scan=11
1250 Factor=False
1260 Interference=False
1270 Numpeak=4
1280 Showdefaults:!
1290 CALL Field_values(X_gdu_max,Y_gdu_max,Numpeak,Interference,Mass(*),Pot(*))
1300 CALL
Parameters(Elem$,Check,Numpeak,Interference,Blockmax,Scan,Factor,Mass(*),Intn(*),Delay(*))
1310 IF Check THEN RETURN
1320 IF NOT Check THEN GOTO Showdefaults
1330 Spikeratios:!
1340 CALL Clear_crt
1350 OUTPUT 1
1360 OUTPUT 1
1370 OUTPUT 1;"1=Ca-Sr Mixed Spike 1 (salt)"
1380 OUTPUT 1;"2=Ca-Sr Mixed Spike 2 (mica)"
1390 OUTPUT 1;"3=Ca43-2(X100)"
1400 OUTPUT 1;"4=Ca43-2"
1410 INPUT "Which spike (type a number then ENTER)?",Sp
1420 IF Sp=1 THEN
1430   Rspike(1)=.013606           !42/43
1440   Rspike(2)=.055219           !42/40
1450   Rspike(3)=.34887           !44/40
1460 END IF
1470 IF Sp=2 THEN
1480   Rspike(1)=.013016
1490   Rspike(2)=.08087
1500   Rspike(3)=.52441
1510 END IF
1520 IF Sp=3 THEN
1530   Rspike(1)=.013135
1540   Rspike(2)=.080413
1550   Rspike(3)=.51197
1560 END IF
1570 IF Sp=4 THEN
1580   Rspike(1)=.013063
1590   Rspike(2)=.082955
1600   Rspike(3)=.53219
1610 END IF
1620 IF Sp<>1 AND Sp<>2 AND Sp<>3 AND Sp<>4 THEN Spikeratios
1630 RETURN
1640 REM Key_handlers
1650 Pause:!
1660 OFF KEY
1670 Halt=True
1680 RETURN
1690 Last_block:!
1700 OFF KEY
1710 Blockmax=Block
1720 RETURN
1730 Escape:!
1740 OFF KEY
1750 CALL Clear_crt
1760 DISP CHR$(131);"PROGRAM ABORTED"

```

```

1770 WAIT 5
1780 GOTO Lastline
1790 New_params: !
1800 OFF KEY
1810 Params=True
1820 RETURN
1830 Summary:!
1840 Index1=2
1850 Index2=5
1860 CALL Final_print(Sample$,Rcorr(*),6,Blockmax)
1870 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Blockmax),4)
1880 CALL Summ_plot(Rcorr(*),Mean(*),Sigma(*),X_gdu_max,Y_gdu_max,Blockmax,Index1,Index2)
1890 DISP CHR$(128)
1900 LOOP
1910   Edit1=0
1920   INPUT "Number of blocks to be edited",Edit1
1930 EXIT IF Edit1=0
1940   CALL Edit(Rcorr(*),Mean(*),Sigma(*),Sigma_int,(Edit1),4,Blockmax)
1950 END LOOP
1960 RETURN
1970 Main:!
1980 OFF KEY
1990 MASS STORAGE IS ":",700,0"
2000 READ LABEL N$ FROM ":",700,0"
2010 A$="DATA"
2020 IF POS(UPC$(N$),A$)=0 THEN
2030   DISP "Insert a data disk in the left drive and press Continue."
2040   PAUSE
2050 END IF
2060 OFF KEY
2061 Today$=DATE$(TIMEDATE)
2062 Now$=TIME$(TIMEDATE)
2063 SELECT Today$[4,6]
2064   CASE "Mar","Jul"
2065     Mon$=Today$[4,4]&Today$[6,6]
2066   CASE ELSE
2067     Mon$=Today$[4,5]
2068 END SELECT
2069 File$=Mon$&Today$[10]&Today$[1,2]&Now$[1,2]
2070 CREATE BDAT File$,52,48
2080 ASSIGN @Data TO File$ !I/O path for disc file
2090 OUTPUT @Data,1;DATE$(TIMEDATE)
2100 OUTPUT @Data,2;Sample$
2110 CALL Start_up(Block$,Sample$,Block,Blockmax,Timbegin)
2120 CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
2130 CALL New_zeros(Baseline(*),Timebase(*),Numpeak,Interference)
2140 WHILE Block<=Blockmax !main program loop
2150   OFF KEY
2160   Halt=False
2170   Params=False
2180   CALL Center_peaks(Qpeak(*),Numpeak,Interference,Mass(*),Pot(*))
2190   Timfirst=INT((TIMEDATE MOD 86400)/60)-Timbegin
2200   Scantime=(SUM(Delay)+SUM(Intn))/60
2210   Timlast=Timfirst+PROUND(Scantime*(Scan+4),0)
2220   CALL Plot(Qpeak(*),X_gdu_max,Y_gdu_max,Timfirst,Timlast)
2230   Block$=VAL$(Block)&" of "&VAL$(Blockmax)
2240   CALL
Measure_peaks(Element$,Elem$,Block$,Baseline(*),Peak(*),Timep(*),Scan,Numpeak,Mass(*),Pot(*),Delay(*),Intn(*),Timbegin,Block,Params)
2250   IF Params THEN GOTO Hold
2260   CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
2270   ON KEY 9 LABEL "ABORT RUN!" GOTO Escape

```

```

2280   ON KEY 5 LABEL "FINISH" GOSUB Last_block
2290   ON KEY 2 LABEL "PAUSE" GOSUB Pause
2300   ON KEY 7 LABEL "Change Params" GOSUB New_params
2310   BEEP 880,2
2320   DISP CHR$(128);"Calculating results"
2330   CALL
Ratios(Traw$(*),Baseline(*),Timebase(*),Timep(*),Peak(*),Netpeak(*),Rraw(*),Avetime(*),Num
peak,Scan,Interference)
2340   CALL
Print_out1(Traw$(*),Baseline(*),Netpeak(*),Rraw(*),Factor,Numpeak,Interference,Block,Scan,
Mass(*),Pot(*))
2350   CALL
Statistics(Avetime(*),Rraw(*),Meanraw(*),Percent_dev(*),Numpeak,Scan,Rejected(*))
2360   CALL Print_out2(Meanraw(*),Percent_dev(*),Numpeak,Rejected(*),Mass(*))
2370   IF Prog=3 THEN
2380     GOSUB Spikesub
2390   ELSE
2400     GOSUB Massdisc
2410   END IF
2420   FOR B=1 TO 6
2430     Rstore(B)=Rcorr(B,Block)
2440   NEXT B
2450   OUTPUT @Data,Block+2;Rstore(*)
2460   CALL Grandstats(Rcorr(*),Mean(*),Sigma(*),Sigma_int,4,(Block))
2470   GOSUB Printout3
2480 Hold:
2490   IF Halt THEN
2500     DISP CHR$(128);"Press CONTINUE to resume data acquisition."
2510     PAUSE
2520   END IF
2530   IF Params THEN
2540     CALL Clear_crt
2550     GOSUB Showdefaults
2560     DISP CHR$(128);"Press CONTINUE to resume data acquisition."
2570     PAUSE
2580   END IF
2590   Block=Block+1
2600   Block$=VAL$(Block)&" of "&VAL$(Blockmax)
2610   IF Params AND Block<=Blockmax THEN CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*
),Intn(*),Pot(*))
2620   CALL New_zeros(Baseline(*),Timebase(*),Numpeak,Interference)
2630 END WHILE
2640 OFF KEY
2650 GOSUB Summary
2660 Lastline: MASS STORAGE IS ":,700,1"
2670 CALL Clear_crt
2680 PRINT USING "@"
2690 LOAD "CONTROL"
2700 END

```

RBBDM

```

10   REM RBBDM                      Brian D. Marshall  7/88
20   OPTION BASE 1
30   PRINTER IS PRT
40   CONTROL 1,4;0                 !Turns off DISPLAY FUNCTIONS
50   LOADSUB ALL FROM "Subfile"
60   DIM
AS[4],Block$[8],Elem$[2],Element$[10],File$[6],N$[6],Sample$[35],Traw$(2,24)[8],Fil$[1]
70   INTEGER Block,Blockmax,Check,Delay(2),Edit1,Factor,False,Halt
80   INTEGER Index1,Index2,Interference,Intn(2),Mass(2),Numpeak,Params,Pot(2)
90   INTEGER Prog,Rejected(2),Scan,Timbegin,Timfirst,Timlast,True
100  REAL Avetime(2,24),Baseline(2,2),Disc,Mean(2)
110  REAL Meanraw(2),Netpeak(2,24),Peak(2,24),Percent_dev(2)
120  REAL Qpeak(2),Rcorr(7,50),Rraw(2,24),Rspike,Rstore(4),Scantime

```

```

130 REAL Sigma(2),Sigma_int,Tcorr(2),Timebase(2,2),Timep(2,24)
140 REAL X_gdu_max,Y_gdu_max
150 CALL Clear_crt
160 False=0
170 True= NOT False
180 Element$="Rubidium"
190 Elem$="RB"
200 Y_gdu_max=100*MAX(1,1/RATIO)
210 X_gdu_max=100*MAX(1,RATIO)
220 CALL Identify(Elem$,Sample$,X_gdu_max,Y_gdu_max)
230 INPUT "Triple filament as perchlorate (T) or single filament as chloride (S)?",Fil$
240 ON KEY 0 LABEL "Sample" GOTO Unspiked
250 ON KEY 2 LABEL "Spiked sample" GOTO Spiked
260 ON KEY 4 LABEL "Spike" GOTO Spike
270 WHILE True
280 END WHILE
290 Unspiked:!
300 Prog=1
310 GOTO Initialize
320 Spiked:!
330 Prog=3
340 GOTO Initialize
350 Spike:!
360 Prog=2
370 Initialize:!
380 OFF KEY
390 GRAPHICS OFF
400 IF Prog=3 THEN GOSUB Spikeratios
410 GOSUB Defaults
420 ON KEY 2 LABEL "START" GOTO Main
430 ON KEY 9 LABEL "Escape" GOTO Initialize
440 ON KEY 5 LABEL "ABORT RUN!" GOTO Escape
450 WHILE True
460 END WHILE
470 Massdisc:!
480 Rcorr(1,Block)=1 !85S/87T
490 Rcorr(2,Block)=Meanraw(2)*(1+2*Disc) !87/85
500 Rcorr(3,Block)=Percent_dev(2)
510 Rcorr(4,Block)=Block
520 Rcorr(7,Block)=1/Rcorr(3,Block)^2
530 RETURN
540 Spikesub:!
550 Tcorr(2)=Meanraw(2)*(1+2*Disc)
560 Rcorr(1,Block)=(1-Tcorr(2)*Rspike)/(Tcorr(2)-.38571)
570 Rcorr(2,Block)=(Tcorr(2)*Rspike-1)/Rcorr(1,Block)+Tcorr(2)
580 Rcorr(3,Block)=Percent_dev(2)
590 Rcorr(4,Block)=Block
600 Rcorr(7,Block)=1/Rcorr(3,Block)^2
610 RETURN
620 Printout3: !
630 Format3: IMAGE 2(4X,3D.7D)
640 IF Prog=3 THEN
650 PRINT
660 PRINT "Spike ratio: ";DROUND(Rspike,6)
670 END IF
680 PRINT
690 PRINT "Mass discrimination corrected (=";Disc;"per a.m.u.)."
700 PRINT
710 PRINT USING "2(9X,7A)";"85S/87T";" 87/85 "
720 PRINT USING Format3;Rcorr(1,Block);Rcorr(2,Block)
730 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Block),2)
740 RETURN
750 Defaults: !
760 Blockmax=10
770 Intn(1)=5
780 Intn(2)=5

```

```

790 Delay(1)=5
800 Delay(2)=5
810 Pot(1)=2040
820 Pot(2)=2735
830 Mass(1)=85
840 Mass(2)=87
850 Scan=11
860 Factor=False
870 Interference=False
880 Numpeak=2
890 IF UPC$(Fil$)="T" THEN
900   Disc=.0017
910 ELSE
920   Disc=.002
930 END IF
940 Showdefaults: !
950 CALL Field_values(X_gdu_max,Y_gdu_max,Numpeak,Interference,Mass(*),Pot(*))
960 CALL
Parameters(Elem$,Check,Numpeak,Interference,Blockmax,Scan,Factor,Mass(*),Intn(*),Delay(*),
Disc)
970 IF Check THEN RETURN
980 IF NOT Check THEN GOTO Showdefaults
990 Spikeratios: !
1000 Rspike=.00833 !85/87
1010 RETURN
1020 REM Key_handlers
1030 Pause:!
1040 OFF KEY
1050 Halt=True
1060 RETURN
1070 Last_block:!
1080 OFF KEY
1090 Blockmax=Block
1100 RETURN
1110 Escape:!
1120 OFF KEY
1130 CALL Clear crt
1140 DISP CHR$(131);"PROGRAM ABORTED"
1150 WAIT 5
1160 GOTO Lastline
1170 New_params:!
1180 OFF KEY
1190 Params=True
1200 RETURN
1210 Summary:!
1220 Index1=(Prog<>3)+1
1230 Index2=3
1240 CALL Final_print(Sample$,Rcorr(*),4,(Blockmax))
1250 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Blockmax),2)
1260 CALL Summ_plot(Rcorr(*),Mean(*),Sigma(*),X_gdu_max,Y_gdu_max,Blockmax,Index1,Index2)
1270 DISP CHR$(128)
1280 LOOP
1290   Edit1=0
1300   INPUT "Number of blocks to be edited",Edit1
1310   EXIT IF Edit1=0
1320   CALL Edit(Rcorr(*),Mean(*),Sigma(*),Sigma_int,(Edit1),2,Blockmax)
1330 END LOOP
1340 RETURN
1350 Main:!
1360 OFF KEY
1370 CALL Start_up(Block$,Sample$,Block,Blockmax,Timbegin)
1380 CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),
Intn(*),Pot(*))
1390 CALL New_zeros(Baseline(*),Timebase(*),Numpeak,Interference)
1400 WHILE Block<=Blockmax

```

```

1410 OFF KEY
1420 Halt=False
1430 Params=False
1440 CALL Center_peaks(Qpeak(*), Numpeak, Interference, Mass(*), Pot(*))
1450 Timfirst=INT((TIMEDATE MOD 86400)/60)-Timbegin
1460 Scantime=(SUM(Delay)+SUM(Intn))/60
1470 Timlast=Timfirst+PROUND(Scantime*(Scan+4), 0)
1480 CALL Plot(Qpeak(*), X_gdu_max, Y_gdu_max, Timfirst, Timlast)
1490 Block$=VAL$(Block)&" of "&VAL$(Blockmax)
1500 CALL
Measure_peaks(Element$, Elem$, Block$, Baseline(*), Peak(*), Timep(*), Scan, Numpeak, Mass(*), Pot(
*), Delay(*), Intn(*), Timbegin, Block, Params)
1510 IF Params THEN GOTO Hold
1520 CALL
Measure_zeros(Block$, Element$, Baseline(*), Timebase(*), Numpeak, Interference, Mass(*), Delay(*
), Intn(*), Pot(*))
1530 ON KEY 9 LABEL "ABORT RUN!" GOTO Escape
1540 ON KEY 5 LABEL "FINISH" GOSUB Last_block
1550 ON KEY 2 LABEL "PAUSE" GOSUB Pause
1560 ON KEY 7 LABEL "Change Params" GOSUB New_params
1570 BEEP 880,2
1580 DISP CHR$(128); "Calculating results"
1590 CALL
Ratios(Traw$(*), Baseline(*), Timebase(*), Timep(*), Peak(*), Netpeak(*), Rraw(*), Avetime(*), Num
peak, Scan, Interference)
1600 CALL
Print_out1(Traw$(*), Baseline(*), Netpeak(*), Rraw(*), Factor, Numpeak, Interference, Block, Scan,
Mass(*), Pot(*))
1610 CALL
Statistics(Avetime(*), Rraw(*), Meanraw(*), Percent_dev(*), Numpeak, Scan, Rejected(*))
1620 CALL Print_out2(Meanraw(*), Percent_dev(*), Numpeak, Rejected(*), Mass(*))
1630 IF Prog=3 THEN
1640 GOSUB Spikesub
1650 ELSE
1660 GOSUB Massdisc
1670 END IF
1680 CALL Grandstats(Rcorr(*), Mean(*), Sigma(*), Sigma_int, 2, (Block))
1690 GOSUB Printout3
1700 Hold:
1710 IF Halt THEN
1720 DISP CHR$(128); "Press CONTINUE to resume data acquisition."
1730 PAUSE
1740 END IF
1750 IF Params THEN
1760 CALL Clear_crt
1770 GOSUB Showdefaults
1780 DISP CHR$(128); "Press CONTINUE to resume data acquisition."
1790 PAUSE
1800 END IF
1810 Block=Block+1
1820 Block$=VAL$(Block)&" of "&VAL$(Blockmax)
1830 IF Params AND Block<=Blockmax THEN CALL
Measure_zeros(Block$, Element$, Baseline(*), Timebase(*), Numpeak, Interference, Mass(*), Delay(*
), Intn(*), Pot(*))
1840 CALL New_zeros(Baseline(*), Timebase(*), Numpeak, Interference)
1850 END WHILE
1860 OFF KEY
1870 GOSUB Summary
1880 Lastline:
1890 CALL Clear_crt
1900 PRINT USING "@"
1910 LOAD "CONTROL"
1920 END

```

SRBDM

10 REM SRBDM

Brian D. Marshall 7/88

```

20  OPTION BASE 1
30  PRINTER IS PRT
40  CONTROL 1,4;0          !Turns off DISPLAY FUNCTIONS.
50  LOADSUB ALL FROM "Subfile"
60  DIM
A$[4],Block$[8],Elem$[2],Element$[10],File$[6],N$[6],Sample$[35],Traw$(4,24)[8],Today$[11],
Now$[8],Mon$[3]
70  INTEGER B,Block,Blockmax,Check,Delay(5),Edit1,Factor,False,Halt,I
80  INTEGER Index1,Index2,Interference,Intn(5),Iter,Mass(5),Numpeak,Params
90  INTEGER Pot(5),Prog,Rejected(4),Scan,Sp,Timbegin,Timfirst,Timlast,True
100 REAL Alpha,Amu(4),Avetime(5,24),Baseline(5,2),Beta,Disc,Ipeak(2),Itime(2)
110 REAL Mean(4),Meancorr(4),Meanraw(4),Netpeak(5,24),P_exp,Peak(5,24)
120 REAL Percent_dev(4),Qpeak(5),Rcorr(7,50),Rraw(5,24),Rspike(3),Rstore(6)
130 REAL Scantime,Sigma(4),Sigma_int,Sum_p,Tcorr(3),Timebase(5,2),Timep(5,24)
140 REAL X_gdu_max,Y_gdu_max
150 CALL Clear_crt
160 OFF KEY
170 DISP CHR$(128)
180 Y_gdu_max=100*MAX(1,1/RATIO)
190 X_gdu_max=100*MAX(1,RATIO)
200 False=0
210 True= NOT False
220 Elem$="SR"
230 Element$="Strontium"
240 CALL Identify(Elem$,Sample$,X_gdu_max,Y_gdu_max)
260 ON KEY 0 LABEL "Sample" GOTO Unspiked
270 ON KEY 2 LABEL "Spiked sample" GOTO Spiked
280 ON KEY 4 LABEL "Spike" GOTO Spike
290 WHILE True
300 END WHILE
310 Unspiked:
320 Prog=1
330 GOTO Initialize
340 Spiked:
350 Prog=3
360 GOTO Initialize
370 Spike:
380 Prog=2
390 Initialize:
400 OFF KEY
410 GRAPHICS OFF
420 IF Prog=3 THEN GOSUB Spikeratios
430 GOSUB Defaults
440 ON KEY 2 LABEL "START" GOTO Main
450 ON KEY 9 LABEL "Escape" GOTO Initialize
460 ON KEY 5 LABEL "ABORT RUN!" GOTO Escape
470 WHILE True
480 END WHILE
490 Massdisc:
500 Sum_p=LOG(.1194/Meanraw(3))/LOG(Amu(3)/Amu(1))
510 IF Prog=2 THEN Sum_p=0          !No correction for spike run
520 Rcorr(1,Block)=1              !86S/84T
530 Rcorr(2,Block)=(Meanraw(2)/Meanraw(3))*(Amu(2)/Amu(3))^Sum_p      !87/86
540 Rcorr(3,Block)=Meanraw(2)*(Amu(2)/Amu(1))^Sum_p                    !87/88
550 Rcorr(4,Block)=(Meanraw(4)/Meanraw(3))*(Amu(4)/Amu(3))^Sum_p      !84/86
560 Rcorr(5,Block)=SQR(2*Percent_dev(3)^2+Percent_dev(2)^2)
570 Rcorr(6,Block)=Block
580 Rcorr(7,Block)=1/Rcorr(5,Block)^2
590 RETURN
600 Spikesub:
610 FOR I=2 TO 4
620 Meancorr(I)=Meanraw(I)
630 NEXT I
640 Sum_p=0
650 Iter=0
660 REPEAT

```

```

670 Alpha=(Meancorr(4)/Meancorr(3)-.056572)/(1-Meancorr(4)/Meancorr(3)*Rspike(1))
680 Beta=1/(1/Meancorr(3)+Alpha*Rspike(1)*(1/Meancorr(3)-1/Rspike(2)))
690 P_exp=LOG(.1194/Beta)/LOG(Amu(3)/Amu(1))
700 Tcorr(1)=1/Alpha
710 Tcorr(2)=(Meancorr(2)/Meancorr(3))*(Amu(2)/Amu(3))^P_exp
720 Tcorr(3)=Meancorr(2)*(Amu(2)/Amu(1))^P_exp
730 Sum_p=Sum_p+P_exp
740 Iter=Iter+1
750 Meancorr(4)=Meancorr(4)*(Amu(4)/Amu(1))^P_exp
760 Meancorr(3)=Tcorr(3)/Tcorr(2)
770 Meancorr(2)=Tcorr(3)
780 UNTIL ABS(P_exp)<1.E-5
790 DISP "Number of iterations was ";Iter
800 Rcorr(1,Block)=Tcorr(1)
810 Rcorr(2,Block)=Tcorr(2)+Alpha*Rspike(1)*(Tcorr(2)-Rspike(3)/Rspike(2))
820 Rcorr(3,Block)=Rcorr(2,Block)*.1194
830 Rcorr(4,Block)=.056572
840 Rcorr(5,Block)=SQRT(2*Percent_dev(3)^2+Percent_dev(4)^2)
850 Rcorr(6,Block)=Block
860 Rcorr(7,Block)=1/Rcorr(5,Block)^2
870 RETURN
880 Printout3: !
890 Disc=((Amu(3)/Amu(1))^Sum_p-1)/2
900 Format3: IMAGE 5(4X,3D.7D)
910 IF Prog=3 THEN
920 PRINT
930 PRINT "Spike ratios:
";DROUND(Rspike(1),6);", ";DROUND(Rspike(2),6);", ";DROUND(Rspike(3),6)
940 PRINT "Number of iterations was ";Iter
950 END IF
960 PRINT
970 IF Interference THEN PRINT "87Rb subtracted; 87Rb/87Sr=
";DROUND(Netpeak(5,Scan)*.387/Netpeak(2,Scan),6)
980 PRINT "Mass discrimination corrected (p=";DROUND(Sum_p,6);",
alpha=";DROUND(Disc,6);")."
990 PRINT
1000 PRINT USING "5(8X,7A)";"86S/84T";" 87/86 ";" 87/88 ";" 84/86 ";"% Error"
1010 PRINT USING
Format3;Rcorr(1,Block);Rcorr(2,Block);Rcorr(3,Block);Rcorr(4,Block);Rcorr(5,Block)
1020 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Block),4)
1030 RETURN
1040 Defaults:!
1050 Blockmax=30
1060 Intn(1)=5
1070 Intn(2)=7
1080 Intn(3)=7
1090 Intn(4)=7
1100 Intn(5)=9
1110 Delay(1)=5
1120 Delay(2)=6
1130 Delay(3)=4
1140 Delay(4)=4
1150 Delay(5)=6
1160 Pot(1)=3025
1170 Pot(2)=2680
1180 Pot(3)=2330
1190 Pot(4)=1630
1200 Pot(5)=1980
1210 Mass(1)=88
1220 Mass(2)=87
1230 Mass(3)=86
1240 Mass(4)=84
1250 Mass(5)=85
1260 Amu(1)=87.905625
1270 Amu(2)=86.908890
1280 Amu(3)=85.909273

```

```

1290 Amu(4)=83.913428
1300 Scan=11
1310 Factor=False
1320 Numpeak=4
1330 Interference=True
1340 Showdefaults:!
1350 CALL Field_values(X_gdu_max,Y_gdu_max,Numpeak,Interference,Mass(*),Pot(*))
1360 CALL
Parameters (Elem$,Check,Numpeak,Interference,Blockmax,Scan,Factor,Mass(*),Intn(*),Delay(*))
1370 IF Check THEN RETURN
1380 IF NOT Check THEN GOTO Showdefaults
1390 Spikeratios:!
1400 CALL Clear_crt
1410 OUTPUT 1
1420 OUTPUT 1
1430 OUTPUT 1;"1=Sr-Ca Mixed Spike 1 (salt)"
1440 OUTPUT 1;"2=Sr-Ca Mixed Spike 2 (mica)"
1450 OUTPUT 1;"3=Sr KF-2"
1460 INPUT "Which spike (type a number then ENTER)?",Sp
1470 IF Sp=1 THEN
1480     Rspike(1)=.003402           !86/84
1490     Rspike(2)=.50446           !86/88
1500     Rspike(3)=.13505           !87/88
1510 END IF
1520 IF Sp=2 THEN
1530     Rspike(1)=.076066
1540     Rspike(2)=5.6316
1550     Rspike(3)=.20517
1560 END IF
1570 IF Sp=3 THEN
1580     Rspike(1)=.0021282
1590     Rspike(2)=.30997
1600     Rspike(3)=.12975
1610 END IF
1620 IF Sp<>1 AND Sp<>2 AND Sp<>3 THEN Spikeratios
1630 RETURN
1640 REM Key_handlers
1650 Pause:!
1660 OFF KEY
1670 Halt=True
1680 RETURN
1690 Last_block:!
1700 OFF KEY
1710 Blockmax=Block
1720 RETURN
1730 Escape:!
1740 OFF KEY
1750 CALL Clear_crt
1760 DISP CHR$(131);"PROGRAM ABORTED"
1770 WAIT 5
1780 GOTO Lastline
1790 New_params:!
1800 OFF KEY
1810 Params=True
1820 RETURN
1830 Summary:!
1840 Index1=2
1850 Index2=5
1860 CALL Final_print(Sample$,Rcorr(*),6,Blockmax)
1870 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Blockmax),4)
1880 CALL Summ_plot(Rcorr(*),Mean(*),Sigma(*),X_gdu_max,Y_gdu_max,Blockmax,Index1,Index2)
1890 DISP CHR$(128)
1900 LOOP
1910     Edit1=0
1920     INPUT "Number of blocks to be edited",Edit1
1930 EXIT IF Edit1=0

```

```

1940 CALL Edit (Rcorr(*), Mean(*), Sigma(*), Sigma_int, (Edit1), 4, Blockmax)
1950 END LOOP
1960 RETURN
1970 Main:
1980 OFF KEY
1990 MASS STORAGE IS ":", 700, 0"
2000 READ LABEL N$ FROM ":", 700, 0"
2010 A$="DATA"
2020 IF POS(UPC$(N$), A$)=0 THEN
2030 DISP "Insert a data disk in the left drive and press Continue."
2040 PAUSE
2050 END IF
2051 Today$=DATE$(TIMEDATE)
2052 Now$=TIME$(TIMEDATE)
2053 SELECT Today$[4, 6]
2054 CASE "Mar", "Jul"
2055 Mon$=Today$[4, 4]&Today$[6, 6]
2056 CASE ELSE
2057 Mon$=Today$[4, 5]
2058 END SELECT
2059 File$=Mon$&Today$[10]&Today$[1, 2]&Now$[1, 2]
2060 CREATE BDAT File$, 52, 48
2070 ASSIGN @Data TO File$ !I/O path for disc file
2080 OUTPUT @Data, 1; DATE$(TIMEDATE)
2090 OUTPUT @Data, 2; Sample$
2100 CALL Start_up(Block$, Sample$, Block, Blockmax, Timbegin)
2110 CALL
Measure_zeros(Block$, Element$, Baseline(*), Timebase(*), Numpeak, Interference, Mass(*), Delay(*), Intn(*), Pot(*))
2120 IF Interference THEN CALL
Measure_inter(Block$, Baseline(*), Ipeak(*), Itime(*), Delay(*), Mass(*), Intn(*), Pot(*))
2130 CALL New_zeros(Baseline(*), Timebase(*), Numpeak, Interference, Ipeak(*), Itime(*))
2140 WHILE Block<=Blockmax !main program loop
2150 OFF KEY
2160 Halt=False
2170 Params=False
2180 CALL Center_peaks(Qpeak(*), Numpeak, Interference, Mass(*), Pot(*))
2190 Timfirst=INT((TIMEDATE MOD 86400)/60)-Timbegin
2200 Scantime=(SUM(Delay)+SUM(Intn))/60-(Delay(5)+Intn(5))/60
2210 Timlast=Timfirst+PROUND(Scantime*(Scan+4), 0)
2220 CALL Plot(Qpeak(*), X_gdu_max, Y_gdu_max, Timfirst, Timlast)
2230 Block$=VAL$(Block)&" of "&VAL$(Blockmax)
2240 CALL
Measure_peaks(Element$, Elem$, Block$, Baseline(*), Peak(*), Timep(*), Scan, Numpeak, Mass(*), Pot(*), Delay(*), Intn(*), Timbegin, Block, Params)
2250 IF Params THEN GOTO Hold
2260 CALL
Measure_zeros(Block$, Element$, Baseline(*), Timebase(*), Numpeak, Interference, Mass(*), Delay(*), Intn(*), Pot(*))
2270 IF Interference THEN CALL
Measure_inter(Block$, Baseline(*), Ipeak(*), Itime(*), Delay(*), Mass(*), Intn(*), Pot(*))
2280 ON KEY 9 LABEL "ABORT RUN!" GOTO Escape
2290 ON KEY 5 LABEL "FINISH" GOSUB Last_block
2300 ON KEY 2 LABEL "PAUSE" GOSUB Pause
2310 ON KEY 7 LABEL "Change Params" GOSUB New_params
2320 BEEP 880, 2
2330 DISP CHR$(128); "Calculating results"
2340 CALL
Ratios(Traw$(*), Baseline(*), Timebase(*), Timep(*), Peak(*), Netpeak(*), Rraw(*), Avetime(*), Numpeak, Scan, Interference, Ipeak(*), Itime(*))
2350 CALL
Print_out1(Traw$(*), Baseline(*), Netpeak(*), Rraw(*), Factor, Numpeak, Interference, Block, Scan, Mass(*), Pot(*))
2360 CALL
Statistics(Avetime(*), Rraw(*), Meanraw(*), Percent_dev(*), Numpeak, Scan, Rejected(*))
2370 CALL Print_out2(Meanraw(*), Percent_dev(*), Numpeak, Rejected(*), Mass(*))

```

```

2380 IF Prog=3 THEN
2390   GOSUB Spikesub
2400 ELSE
2410   GOSUB Massdisc
2420 END IF
2430 FOR B=1 TO 6
2440   Rstore(B)=Rcorr(B,Block)
2450 NEXT B
2460 OUTPUT @Data,Block+2;Rstore(*)
2470 CALL Grandstats(Rcorr(*),Mean(*),Sigma(*),Sigma_int,4,(Block))
2480 GOSUB Printout3
2490 Hold:
2500 IF Halt THEN
2510   DISP CHR$(128);"Press CONTINUE to resume data acquisition."
2520   PAUSE
2530 END IF
2540 IF Params THEN
2550   CALL Clear_crt
2560   GOSUB Showdefaults
2570   DISP CHR$(128);"Press CONTINUE to resume data acquisition."
2580   PAUSE
2590 END IF
2600 Block=Block+1
2610 Block$=VAL$(Block)&" of "&VAL$(Blockmax)
2620 IF Params AND Block<=Blockmax THEN
2630   CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
2640   IF Interference THEN CALL
Measure_inter(Block$,Baseline(*),Ipeak(*),Itime(*),Delay(*),Mass(*),Intn(*),Pot(*))
2650   END IF
2660   CALL New_zeros(Baseline(*),Timebase(*),Numpeak,Interference,Ipeak(*),Itime(*))
2670 END WHILE
2680 OFF KEY
2690 GOSUB Summary
2700 Lastline: MASS STORAGE IS ":,700,1"
2710 CALL Clear_crt
2720 PRINT USING "@ "
2730 LOAD "CONTROL"
2740 END

```

IRBDM

```

10 REM IRBDM Brian D. Marshall 7/88
20 OPTION BASE 1
30 PRINTER IS CRT
40 CONTROL 1,4;0 !Turns off DISPLAY FUNCTIONS
50 LOADSUB ALL FROM "Subfile"
60 DIM A$(4),Block$(8),Elem$(2),Element$(10),File$(6),N$(6),Sample$(35),Traw$(5,24)[8]
70 INTEGER Block,Blockmax,Check,Delay(5),Edit1,Factor,False,Halt
80 INTEGER Index1,Index2,Interference,Intn(5),Iter,Mass(5),Numpeak,Params
90 INTEGER Pot(5),Rejected(5),Scan,Timbegin,Timfirst,Timlast,True
100 INTEGER P_mass(5),P_pot(5),Sel,Numsel
110 REAL Avetime(5,24),Baseline(5,2),Mean(5)
120 REAL Meanraw(5),Netpeak(5,24),Peak(5,24),Percent_dev(5)
130 REAL Qpeak(5),Rcorr(7,50),Rraw(5,24),Scantime
140 REAL Sigma(5),Sigma_int,Timebase(5,2),Timep(5,24)
150 REAL X_gdu_max,Y_gdu_max
160 CALL Clear_crt
170 Y_gdu_max=100*MAX(1,1/RATIO)
180 X_gdu_max=100*MAX(1,RATIO)
190 False=0
200 True= NOT False
210 Elem$="IR"
220 CALL Identify(Elem$,Sample$,X_gdu_max,Y_gdu_max)
230 Element:

```

```
240 OFF KEY
250 INPUT "Element to be analyzed (U, TH, PB, RE, RB, K, SR, CA)?",El$
260 Elem$=UPC$(El$)
270 SELECT Elem$
280 CASE ="U"
290   Element$="Uranium"
300   P_mass(1)=233
310   P_mass(2)=234
320   P_mass(3)=235
330   P_mass(4)=236
340   P_mass(5)=238
350   P_pot(1)=1000
360   P_pot(2)=1500
370   P_pot(3)=2000
380   P_pot(4)=2500
390   P_pot(5)=3000
400   Numsel=5
410   GOSUB Select
420 CASE ="TH"
430   Element$="Thorium"
440   Numpeak=2
450   REDIM Intn(2),Delay(2),Mass(2),Pot(2)
460   Mass(1)=230
470   Mass(2)=232
480   Pot(1)=700
490   Pot(2)=1350
500 CASE ="RB"
510   Element$="Rubidium"
520   Numpeak=2
530   REDIM Intn(2),Delay(2),Mass(2),Pot(2)
540   Mass(1)=85
550   Mass(2)=87
560   Pot(1)=1952
570   Pot(2)=2655
580 CASE ="K"
590   Element$="Potassium"
600   P_mass(1)=39
610   P_mass(2)=40
620   P_mass(3)=41
630   P_pot(1)=900
640   P_pot(2)=1350
650   P_pot(3)=1860
660   Numsel=3
670   GOSUB Select
680 CASE ="RE"
690   Element$="Rhenium"
700   Numpeak=2
710   REDIM Intn(2),Delay(2),Mass(2),Pot(2)
720   Mass(1)=185
730   Mass(2)=187
740   Pot(1)=1000
750   Pot(2)=2000
760 CASE ="PB"
770   Element$="Lead"
780   P_mass(1)=204
790   P_mass(2)=205
800   P_mass(3)=206
810   P_mass(4)=207
820   P_mass(5)=208
830   P_pot(1)=1000
840   P_pot(2)=1500
850   P_pot(3)=2000
860   P_pot(4)=2500
870   P_pot(5)=3000
880   Numsel=5
890   GOSUB Select
```

```

900 CASE "SR"
910   Element$="Strontium"
920   P_mass(1)=84
930   P_mass(2)=86
940   P_mass(3)=87
950   P_mass(4)=88
960   P_pot(1)=1545
970   P_pot(2)=2260
980   P_pot(3)=2610
990   P_pot(4)=2960
1000  Numsel=4
1010  GOSUB Select
1020 CASE "CA"
1030   Element$="Calcium"
1040   P_mass(1)=40
1050   P_mass(2)=42
1060   P_mass(3)=43
1070   P_mass(4)=44
1080   P_mass(5)=48
1090   P_pot(1)=1700
1100   P_pot(2)=2600
1110   P_pot(3)=3040
1120   P_pot(4)=3475
1130   P_pot(5)=4000
1140   Numsel=5
1150   GOSUB Select
1160 CASE ELSE
1170   GOTO Element
1180 END SELECT
1190 CALL Clear_crt
1200 PRINT
1210 PRINT "Selected ratios:"
1220 FOR R=2 TO Numpeak
1230   PRINT
1240   PRINT Mass(R);"/";Mass(1)
1250 NEXT R
1260 ON KEY 2 LABEL "OK" GOTO Initialize
1270 ON KEY 9 LABEL "Re-select" GOTO Element
1280 WHILE True
1290 END WHILE
1300 Select:
1310 CALL Clear_crt
1320 INPUT "Number of isotopes to be measured?",Numpeak
1330 IF Numpeak<2 OR Numpeak>5 THEN GOTO Element
1340 REDIM Intn(Numpeak),Delay(Numpeak),Pot(Numpeak),Mass(Numpeak)
1350 PRINT
1360 PRINT
1370 FOR I=1 TO Numsel
1380   PRINT I;"=";P_mass(I)
1390 NEXT I
1400 INPUT "Select reference isotope (type a number then ENTER)?",Sel
1410 IF Sel<1 OR Sel>Numsel THEN GOTO Element
1420 PRINT TABXY(10,Sel+2);"Peak 1 (reference)";
1430 Mass(1)=P_mass(Sel)
1440 Pot(1)=P_pot(Sel)
1450 FOR N=2 TO Numpeak
1460   INPUT "Select next isotope to be measured (type a number then ENTER)?",Sel
1470   IF Sel<1 OR Sel>Numsel THEN GOTO Element
1480   PRINT TABXY(10,Sel+2);"Peak ";N;
1490   Mass(N)=P_mass(Sel)
1500   Pot(N)=P_pot(Sel)
1510 NEXT N
1520 RETURN
1530 Initialize:
1540 PRINTER IS PRT
1550 OFF KEY

```

```

1560 GRAPHICS OFF
1570 GOSUB Defaults
1580 ON KEY 2 LABEL "START" GOTO Main
1590 ON KEY 9 LABEL "Escape" GOTO Initialize
1600 ON KEY 5 LABEL "ABORT RUN!" GOTO Escape
1610 WHILE True
1620 END WHILE
1630 Massdisc:~
1640 Rcorr(1,Block)=Meanraw(2)
1650 Rcorr(2,Block)=Meanraw(3)
1660 Rcorr(3,Block)=Meanraw(4)
1670 Rcorr(4,Block)=Meanraw(5)
1680 Rcorr(5,Block)=Percent_dev(2)
1690 Rcorr(6,Block)=Block
1700 Rcorr(7,Block)=1/Rcorr(5,Block)^2
1710 RETURN
1720 Printout3:~
1730 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Block),Numpeak-1)
1740 RETURN
1750 Defaults:~
1760 Blockmax=10
1770 FOR N=1 TO Numpeak
1780   Intn(N)=5
1790   Delay(N)=5
1800 NEXT N
1810 Scan=11
1820 Factor=False
1830 Interference=False
1840 Showdefaults:~
1850 CALL Field_values(X_gdu_max,Y_gdu_max,Numpeak,Interference,Mass(*),Pot(*))
1860 CALL
Parameters(Elem$,Check,Numpeak,Interference,Blockmax,Scan,Factor,Mass(*),Intn(*),Delay(*))
1870 IF Check THEN RETURN
1880 IF NOT Check THEN GOTO Showdefaults
1890 REM Key_handlers
1900 Pause:~
1910 OFF KEY
1920 Halt=True
1930 RETURN
1940 Last_block:~
1950 OFF KEY
1960 Blockmax=Block
1970 RETURN
1980 Escape:~
1990 OFF KEY
2000 CALL Clear crt
2010 DISP CHR$(131);"PROGRAM ABORTED"
2020 WAIT 5
2030 GOTO Lastline
2040 New_params:~
2050 OFF KEY
2060 Params=True
2070 RETURN
2080 Summary:~
2090 Index1=1
2100 Index2=5
2110 CALL Final_print(Sample$,Rcorr(*),6,Blockmax)
2120 CALL Print_means(Mean(*),Sigma(*),Sigma_int,(Blockmax),Numpeak-1)
2130 CALL Summ_plot(Rcorr(*),Mean(*),Sigma(*),X_gdu_max,Y_gdu_max,Blockmax,Index1,Index2)
2140 DISP CHR$(128)
2150 LOOP
2160   Edit1=0
2170   INPUT "Number of blocks to be edited",Edit1
2180 EXIT IF Edit1=0
2190   CALL Edit(Rcorr(*),Mean(*),Sigma(*),Sigma_int,(Edit1),Numpeak-1,Blockmax)
2200 END LOOP

```

```

2210 RETURN
2220 Main:
2230 OFF KEY
2240 PRINTER IS PRT
2250 CALL Start_up(Block$,Sample$,Block,Blockmax,Timbegin)
2260 CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
2270 CALL New_zeros(Baseline(*),Timebase(*),Numpeak,Interference)
2280 WHILE Block<=Blockmax
2290   OFF KEY
2300   Halt=False
2310   Params=False
2320   CALL Center_peaks(Qpeak(*),Numpeak,Interference,Mass(*),Pot(*))
2330   Timfirst=INT((TIMEDATE MOD 86400)/60)-Timbegin
2340   Scantime=(SUM(Delay)+SUM(Intn))/60
2350   Timlast=Timfirst+PROUND(Scantime*(Scan+4),0)
2360   CALL Plot(Qpeak(*),X_gdu_max,Y_gdu_max,Timfirst,Timlast)
2370   Block$=VAL$(Block)&" of "&VAL$(Blockmax)
2380   CALL
Measure_peaks(Element$,Elem$,Block$,Baseline(*),Peak(*),Timep(*),Scan,Numpeak,Mass(*),Pot(*),Delay(*),Intn(*),Timbegin,Block,Params)
2390   IF Params THEN GOTO Hold
2400   CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
2410   ON KEY 9 LABEL "ABORT RUN!" GOTO Escape
2420   ON KEY 5 LABEL "FINISH" GOSUB Last_block
2430   ON KEY 2 LABEL "PAUSE" GOSUB Pause
2440   ON KEY 7 LABEL "Change Params" GOSUB New_params
2450   BEEP 880,2
2460   DISP CHR$(128);"Calculating results"
2470   CALL
Ratios(Traw$(*),Baseline(*),Timebase(*),Timep(*),Peak(*),Netpeak(*),Rraw(*),Avetime(*),Numpeak,Scan,Interference)
2480   CALL
Print_out1(Traw$(*),Baseline(*),Netpeak(*),Rraw(*),Factor,Numpeak,Interference,Block,Scan,Mass(*),Pot(*))
2490   CALL
Statistics(Avetime(*),Rraw(*),Meanraw(*),Percent_dev(*),Numpeak,Scan,Rejected(*))
2500   CALL Print_out2(Meanraw(*),Percent_dev(*),Numpeak,Rejected(*),Mass(*))
2510   GOSUB Massdisc
2520   CALL Grandstats(Rcorr(*),Mean(*),Sigma(*),Sigma_int,Numpeak-1,(Block))
2530   GOSUB Printout3
2540 Hold:
2550   IF Halt THEN
2560     DISP CHR$(128);"Press CONTINUE to resume data acquisition."
2570     PAUSE
2580   END IF
2590   IF Params THEN
2600     CALL Clear_crt
2610     GOSUB Showdefaults
2620     DISP CHR$(128);"Press CONTINUE to resume data acquisition."
2630     PAUSE
2640   END IF
2650   Block=Block+1
2660   Block$=VAL$(Block)&" of "&VAL$(Blockmax)
2670   IF Params AND Block<=Blockmax THEN CALL
Measure_zeros(Block$,Element$,Baseline(*),Timebase(*),Numpeak,Interference,Mass(*),Delay(*),Intn(*),Pot(*))
2680   CALL New_zeros(Baseline(*),Timebase(*),Numpeak,Interference)
2690 END WHILE
2700 OFF KEY
2710 GOSUB Summary
2720 Lastline:
2730 CALL Clear_crt

```

```
2740 PRINT USING "@"  
2750 LOAD "CONTROL"  
2760 END
```

References

- Dodson M.H. (1978) A linear method for second-degree interpolation in cyclical data collection. *J. Phys. E: Sci. Instrum.* **11**, 296.
- Taylor J.R. (1982) *An introduction to error analysis: the study of uncertainties in physical measurements*. University Science Books, Mill Valley, 270 p.
- Wasserburg G.J., Jacobsen S.B., DePaolo D.J., McCulloch M.T. and Wen T. (1981) Precise determination of Sm/Nd ratios, Sm and Nd isotopic abundances in standard solutions. *Geochim. Cosmochim. Acta* **45**, 2311-2323.